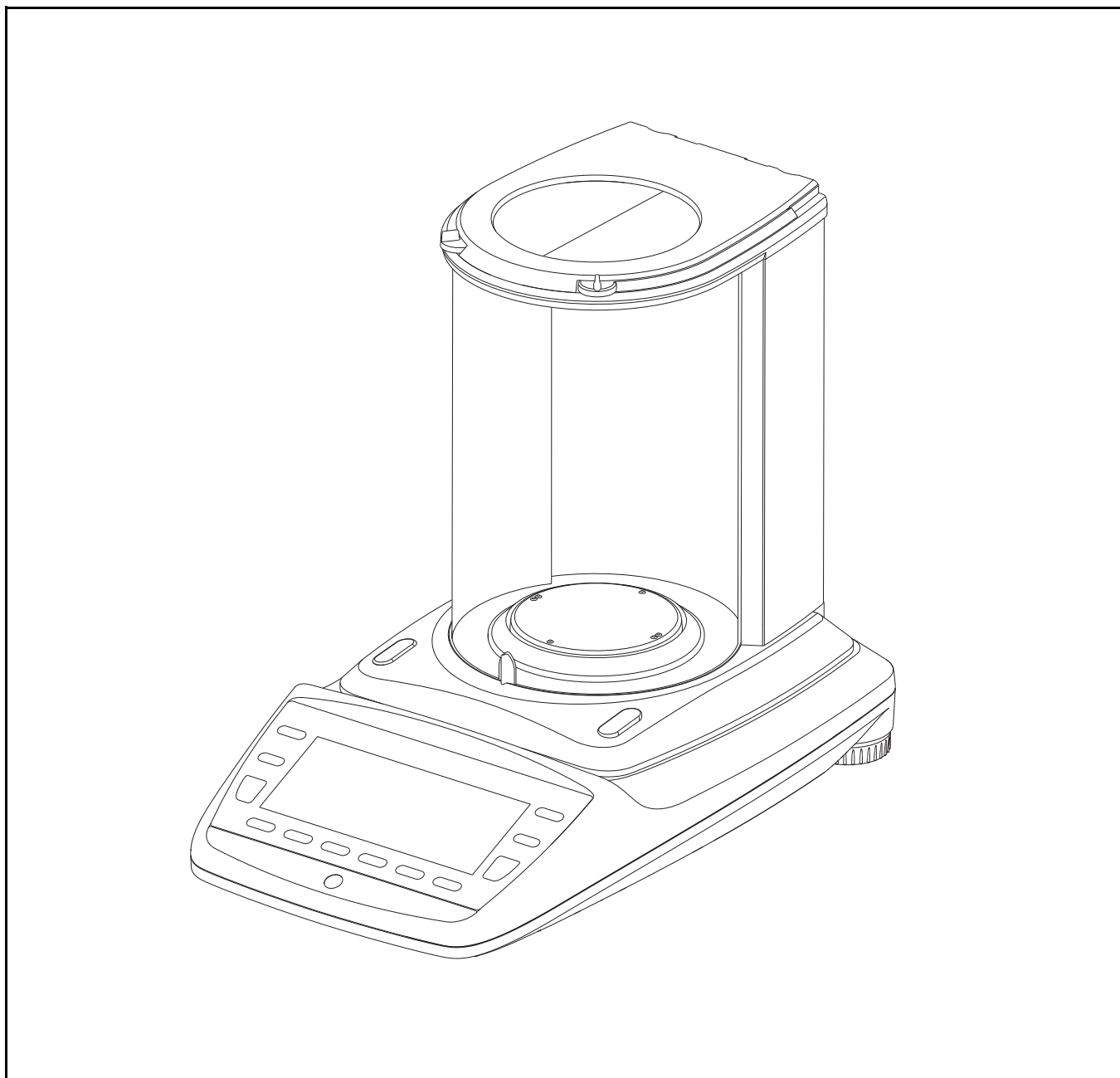
# Precisa 360 EP

Brief Operating Instructions  
Kurzbedienungsanleitung  
Mode d'emploi court  
Manual de Instrucciones corta

## Declaration of conformity

**Declaration of conformity for apparatus with CE mark**  
**Konformitätserklärung für Geräte mit CE-Zeichen**  
**Déclaration de conformité pour appareils portant la marque CE**  
**Declaración de conformidad para aparatos con disitintivo CE**  
**Dichiarazione di conformità per apparecchi contrassegnati con la marcatura CE**

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**Electronic Balance: Precisa Series 360 EP**  
**Precisa Series 360 ES**  
**Precisa Series 321 LS**  
**Precisa Series 321 LX**  
**Precisa Series 321 LT**

EU Directive	Standards	EC-type-approval	Issued by
2014/31/EU	EN 45501: 2015	D11-09-010 T7911	PTB NMI

Date: 20.04.2016

Signature:

R. Grolimund R & D Manager

Precisa Gravimetrics AG , Moosmattstrasse 32 , Postfach 352 , CH-8953 Dietikon

# 1 Safety Instructions and Conformity

## 1.1 Conformity

The device has been manufactured and tested in accordance with the standards and directives set out in the enclosed certificate of conformity.

## 1.2 Safety Instructions

- When using the device in surroundings with increased safety requirements, the corresponding regulations must be observed.
- The device may only be operated with the included power adaptor or power cable.
- Be sure to hold the plug, not the cable, when disconnecting the monitor from an electrical outlet.
- If the power cable, the power adaptor or its cable is damaged, unplug the device immediately from the electrical outlet and keep it from inadvertent operation. The device may only be operated with a power adaptor or a power cable in perfect conditions.
- If there is any reason to believe that it is no longer possible to operate the device without danger, it must be immediately unplugged from the electrical outlet and kept from inadvertent operation.
- In carrying out maintenance work, it is essential to heed the instructions in the chapter "Maintenance and servicing" of the operating instructions.
- Do not operate the device in an area subject to explosion risks or caustic atmospheres.
- Take care when handling with liquids, that no liquid is split into the inside of the balance, on connections of the device, on the power cable or the power adaptor. Should this although happen, unplug the device immediately from the electrical outlet and keep it from inadvertent operation. The device may only be operated again after a re-check made by a Precisa service technician.
- The device may only be used for weighing of solid materials, of liquids filled into secure containers, and for running applications designed especially for the device (e.g. animal weighing or density determinations).
- Do not exceed the maximum permissible load of the balance.
- When using the device in combination with other appliances made by Precisa as well as with appliances produced by other manufacturers, the current regulations for the safe use of the relevant attachments and their application in accordance with instructions must always be observed.

### **DANGER**

**Before connecting the device to the mains, be sure that the operating voltage stated on the device and on the power adaptor or power cable agrees with the local mains voltage. If it doesn't, the equipment must not be connected to the mains at all! In this case contact the Precisa Customer Service.**

### **NOTE**

**If any damage or injury occurs, liability and responsibility rest with the user.**

## 2 Discover the Capabilities of the New Series 360 EP

Work through the following example in order to explore some of the many new features of your new Precisa balance of the series 360 EP. Follow the instructions and keep observing the display.

### Example:

As part of a modification kit, destined to 15 of your customers, you need to pack screws. Each kit is to consist of 126 screws. Since the screws are small and easy to get lost, you round the number up to 130 pieces, in order to keep possible trouble away from your customers. It although doesn't matter, if somebody gets one or two screws more or less. Let, for a better convenience, the balance do the counting of the weighed packs. After all, for your statistics, you would like to know, how many screws you have packed and how accurate has been your working.

### 2.1 1<sup>st</sup> Step

Switch-on the balance:

- Press «**ON/OFF**»

During the start-up sequence the balance executes a self check routine and stays, after about five seconds, ready in the same application before its last switching-off.

In order to follow this example exactly, you may set the factory settings to the balance, but consider, that eventually made previous settings will be lost, wherefore a question window will ask you to confirm:

- Hold «**MENU**»
- Keep pressing {→} until „Settings“ in the headline is selected
- Press {▲} beneath the Administrator icon
- Keep pressing {↓} until the menu item „Set factory settings“ is selected and
- Press {↵} to confirm your selection
- Affirm the question „Set factory settings?“ by pressing {Yes}

Now select your language:

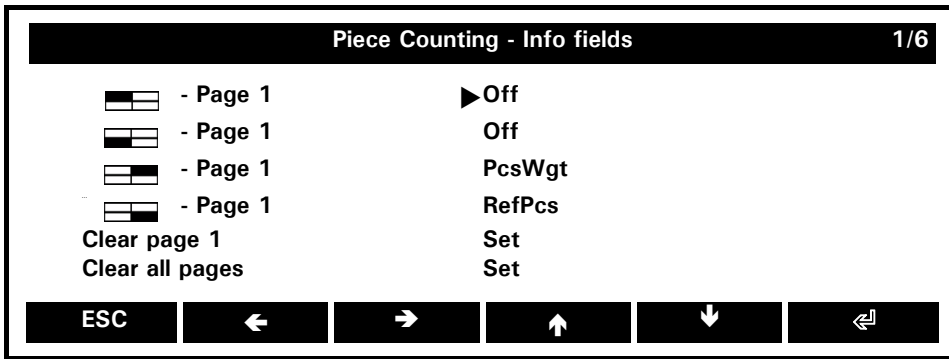
- Hold «**MENU**»
- Keep pressing {→} until „Settings“ in the headline is selected
- Press {▲} beneath the Configuration icon
- Keep pressing {↓} until the menu item „Device settings“ is selected
- Press {↵} to confirm your selection
- Keep pressing {↓} until the menu item „Language“ is selected
- Press {↵} to confirm your selection
- Keep pressing {↓} until in the list box your language is selected and
- Press {↵} to confirm your selection

Return to the application „Weighing“:

- Hold «**MENU**»
- Press {▲} beneath the Weighing icon


Now switch the menu item „Info fields“:

As the CheckWeigher is already enabled, some basic additional information, such as PcsWgt and RefPcs, is going to be displayed in the info page. In our example we want to take a closer look over the shoulder of the balance and let it show us the last and the penultimate stored value too.

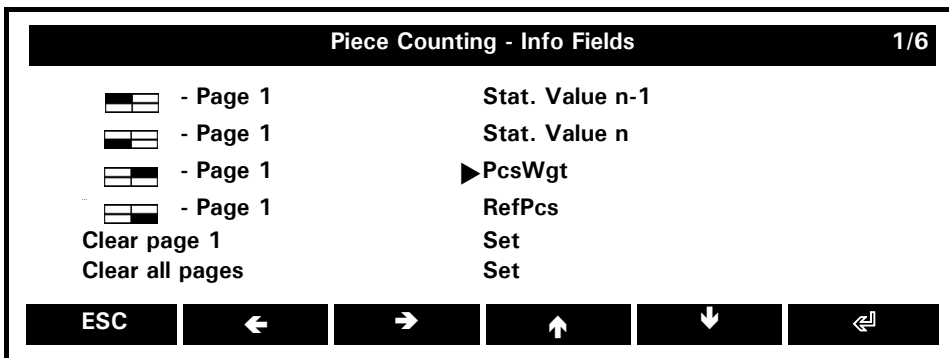Assign the first two info page positions:

English

Return to the application Piece Counting:

- Press {ESC} twice

## ■ 2 Discover the Capabilities of the New Series 360 EP

### 2.3 3<sup>rd</sup> Step

Put the container in which you want to count the screws on the weighing pan and tare the balance:

- Press «0/T»

At first, put exactly 5 screws into the container.

- Press {5 PCS}

The balance takes the quantity showing it in the primary display, while the weight is visible in the secondary display. There appears a rotate-symbol  on its right end which indicates the possibility of a content interchange between the primary and the secondary display by pressing «ROTATE».

In the Info Area the balance is showing under "PcsWgt =" the calculated weight of one screw. Each time the number of pieces increases for at least three, the balance, being stable, re-calculates the reference weight of one piece. During this procedure appears a circle "O" on the left end of the primary display. As soon as it disappears, the new, just calculated piece weight is displayed in its info field. In order to allow the balance to count error-free, it is, at every first weighing-in, advisable to pause for some seconds each time the number of pieces doubles the last calculated one (thus in our case at 10, 20, 40 and 80).

As soon as the first lot of 130 pcs., with the graphic and acoustic aid of the CheckWeigher, is weighed-in, you may store it:

- Press {STO 1}

The function "STO 1" changes to "STO 2", "STO 3" (observe both info fields on the left side of the info area). It stores and counts every weighing-in. Once the last weighing-in is stored, in our case the 15th, it changes to "ENDED 15".

All statistic data are continuously refreshed and available during the entire procedure, as well individually displayed in the info area (remember the hold «ROTATE» -key for running over the info pages), as recallable with {Show Stat.} in an info window.

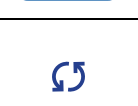
By pressing {Clr Stat.} you may anytime clear all statistic data and restart your weighing-in with the existing procedure settings.

## 3 Operation

### 3.1 The Keys

#### 3.1.1 Functions of the Keys if Operated in an Application

Key	Designation	Function in an Application
	«ON/OFF»	<ul style="list-style-type: none"> <li>• Turns the balance on and off.</li> </ul>
	«PRINT»	<ul style="list-style-type: none"> <li>• Transmits the current measuring result or the content of an info window via interface to the peripherals.</li> </ul>
	«MENU»	<ul style="list-style-type: none"> <li>• Pressing leads to the context menu of the currently executed application.</li> <li>• Holding leads to the home screen from any point.</li> </ul>
	«ROTATE»	<ul style="list-style-type: none"> <li>• Pressing (only if adverted by the rotate symbol  in the activity area) exchanges the measuring results of the primary and the secondary display.</li> <li>• Holding calls up the next info page, continuously holding scrolls cyclically all info pages.</li> </ul>
	«O/T»	<ul style="list-style-type: none"> <li>• Pressing re-zeroes, respectively tares the balance.</li> <li>• Holding leads to the calibration menu.</li> </ul>
	{SOFT KEY}	<ul style="list-style-type: none"> <li>• Each of the six soft keys executes its assigned function, displayed in the function bar, straight above the key itself. If displayed, pressing {→} or {←} calls up further softkey-functions.</li> </ul>

#### 3.1.2 Functions of the Keys if Operated in a Menu

Key	Designation	Function in a Menu
	«ON/OFF»	<ul style="list-style-type: none"> <li>• Turns the balance on and off.</li> </ul>
	«PRINT»	<ul style="list-style-type: none"> <li>• Not functional.</li> </ul>
	«MENU»	<ul style="list-style-type: none"> <li>• Pressing leaves a menu.</li> <li>• Holding leads to the home screen from any point.</li> </ul>
	«ROTATE»	<ul style="list-style-type: none"> <li>• Pressing (only if adverted by the rotate symbol  in the activity area) calls up further soft key functions.</li> </ul>
	«O/T»	<ul style="list-style-type: none"> <li>• Not functional.</li> </ul>

## ■ 3 Operation

Key	Designation	Function in a Menu
	{SOFT KEY}	<ul style="list-style-type: none"> <li>Each of the six soft keys executes its assigned function, displayed in the function bar, straight above the key itself. If available (displayed rotate symbol  in the activity area), pressing «<b>ROTATE</b>» calls up further soft key functions.</li> </ul>

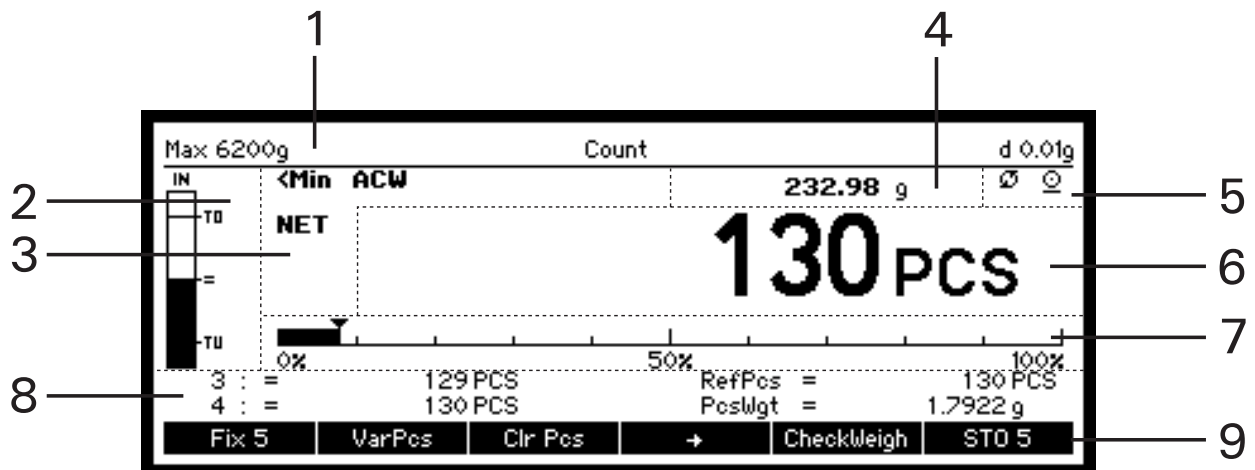
### 3.1.3 Soft Key Functions for Navigation and Data Entry

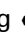
{←}	{left}	Moves the cursor left / scrolls pages
{→}	{right}	Moves the cursor right / scrolls pages
{↑}	{up}	Moves the cursor up / scrolls up lines
{↓}	{down}	Moves the cursor down / scrolls down lines
{▲}	{switch}	Switches the icon-figured function above
{↵}	{Enter}	Selects a menu item / finishes and saves an input
{ESC}	{Escape}	Quits an input without saving the changes
{ ←}	{Backspace}	Deletes the character located on the left side of the cursor („_“)
{DEL}	{Delete}	Deletes the character located above the cursor („_“)
{ABCDEFG}	{Symbol entry}	Keep pressing until the desired symbol appears, e.g. twice for „B”. The same procedure is valid for all symbol entry functions, e.g. {+ -* / = & %} or {67890}
{_}	{Space}	Interword separation

### 3.1.4 Soft Keys and Soft Key Assignments

A soft key is a soft-coded key, a button, located alongside the display, which performs a function displayed near or right above it in the function bar at that moment. (In opposition to that, a hard key is a hard-coded key, which performs just its firmly assigned function, e.g. «**ON/OFF**».) In each application you may assign a choice of various functions to the soft keys. If you assign more than 6 functions, a function {←} or {→} is displayed to switch from the current selection to the next.

### 3.2 The Display (as it appears during a running application)


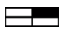
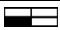
No.	Name of Display Section	Description
1	Headline	<ul style="list-style-type: none"> <li>Names the running application with the weighing range and readability, or the running application and the system clock with date and time.</li> </ul>
2	CheckWeigher	<ul style="list-style-type: none"> <li>This smart aid supports you graphically, and by wish also acoustically, while weighing in.</li> </ul>
3	Notice Area	<ul style="list-style-type: none"> <li>Displays additional information about the measuring result, e.g. „NET“.</li> <li>Advises by displaying „&lt;Min“ of a minimal weight underrun.</li> <li>Advises by displaying „ACW“ (Automatic calibration warning) the balance needs to be calibrated.</li> </ul>
4	Secondary Display	<ul style="list-style-type: none"> <li>Displays the measuring result, if chosen, in another unit, e.g. „PCS“. Observe that the unit symbol only appears at the time the measuring result is stable.</li> </ul>
5	Activity Area	<ul style="list-style-type: none"> <li>A displayed rotate symbol  adverts that by pressing «RO-TATE» the measuring results of the primary and the secondary display can be exchanged.</li> <li>A displayed print symbol adverts that the balance is busy transmitting the current measuring result or the content of an info window via interface to the peripherals.</li> </ul>
6	Primary Display	<ul style="list-style-type: none"> <li>Displays the measuring result in the chosen basic unit. Observe that the unit symbol only appears at the time the measuring result is stable.</li> </ul>
7	Capacity Indicator	<ul style="list-style-type: none"> <li>Displays graphically the rate of the currently utilised weighing range „0% ... 50% ... 100%“.</li> </ul>
8	Info Page Area	<ul style="list-style-type: none"> <li>Displays the selected info page, consisting of up to four assigned info fields available in the menu of the running application. This is to show information like tolerance limits when you are weighing in or various statistical data.</li> </ul>
9	Function Bar	<ul style="list-style-type: none"> <li>Displays straight above every single soft key its, at the moment assigned, function.</li> </ul>

## ■ 3 Operation
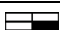
### 3.2.1 Info Pages and Info Fields

There is a screen section called „Info Page Area“ to display application specific information. Each info page has 4 info fields: top left, bottom left, top right and bottom right. In an application, holding the «**ROTATE**» key calls up the next info page and continuously holding it scrolls cyclically all info pages.

**Available info fields (quadrants) for displaying information on the info pages:**

 Page 1 / Top left	 Page 1 / Top right
 Page 1 / Bottom left	 Page 1 / Bottom right
 Page 2 / Top left	 Page 2 / Top right
 Page 2 / Bottom left	 Page 2 / Bottom right
 Page 3 / Top left	 Page 3 / Top right
 Page 3 / Bottom left	 Page 3 / Bottom right
...	

Some applications already display the most usual information, but you may reconfigure the info pages at any time. Therefore switch the menu item Info fields in the context menu of the application and assign the info fields individually.

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