
SWISS MADE 

Series 321

Precisa Gravimetrics AG

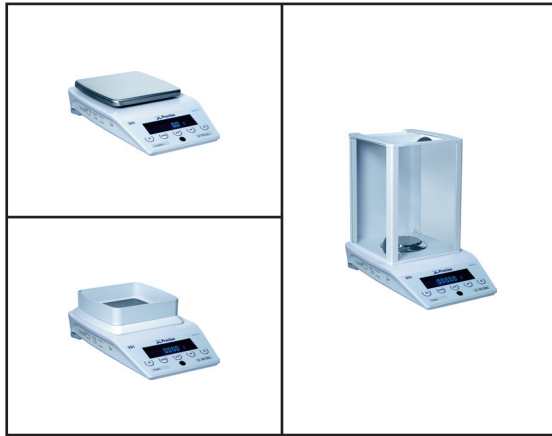Brief instructions / Kurzanleitung / Courte notice d'utilisation / Instrucciones breves


Precisa Gravimetrics AG
Moosmattstrasse 32
CH-8953 Dietikon
Tel.: +41-44-744 28 28
Fax.: +41-44-744 28 38
Email: service@precisa.ch

<http://www.precisa.com>

Customer service

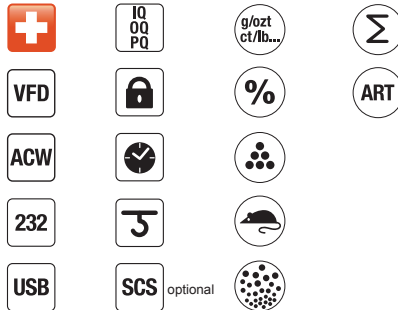Information on and addresses for local customer service centres can be found on our website.

Copyright

These operating instructions are protected by copyright. All rights reserved. No part of these operating instructions may be reproduced, edited, copied or distributed in any form either through photocopying, microfiche, reprinting or any other method, and in particular via any electronic means, without written permission.

© Precisa Gravimetrics AG, 8953 Dietikon, Switzerland, 2013.

Table of contents

| | | | | | |
|-----------|----------------------------------|-----------|------------|---|-----------|
| 1. | Introduction | 6 | 7. | Key menus | 16 |
| 2. | Safety precautions | 6 | 7.1 | Tare menu | 16 |
| 3. | Technical data | 7 | 7.1.1 | Taring the balance..... | 16 |
| 4. | Starting up | 8 | 7.1.2 | Other functions | 16 |
| 4.1 | Choice of a suitable location .. | 8 | 7.2 | Mode menu | 17 |
| 4.2 | Unpacking the balance | 8 | 7.2.1 | Switch to an activated application or back to weighing mode..... | 17 |
| 4.3 | Assembling the balance..... | 9 | 7.2.2 | Other functions | 17 |
| 4.4 | Under-floor weighing | 10 | 7.3 | Print menu | 18 |
| 4.5 | Levelling the balance..... | 11 | 7.3.1 | Print individual value / report..... | 18 |
| 4.6 | Calibrating the balance | 11 | 7.3.2 | Other functions | 18 |
| 5. | Key functions..... | 12 | | | |
| 6. | Overview of menus | 13 | | | |
| 6.1 | Menu control | 13 | | | |
| 6.2 | Configuration menu..... | 14 | | | |
| 6.3 | Application menu..... | 15 | | | |

**DANGER**

Warning of a possible danger which can lead to death or serious physical injuries.

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**WARNING**

Warning of a possibly dangerous situation which can lead to less severe injuries or damage.

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**NOTE**

Tips and important rules on the correct operation of the balance.

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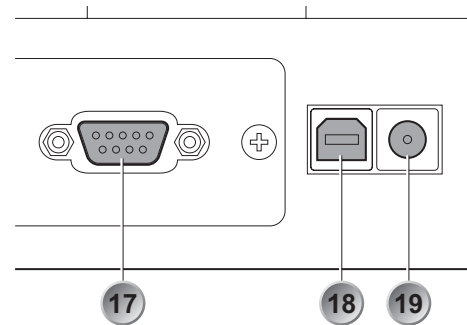
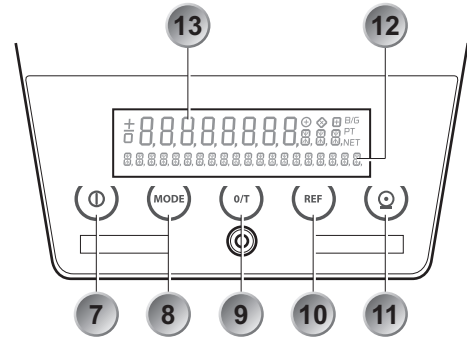
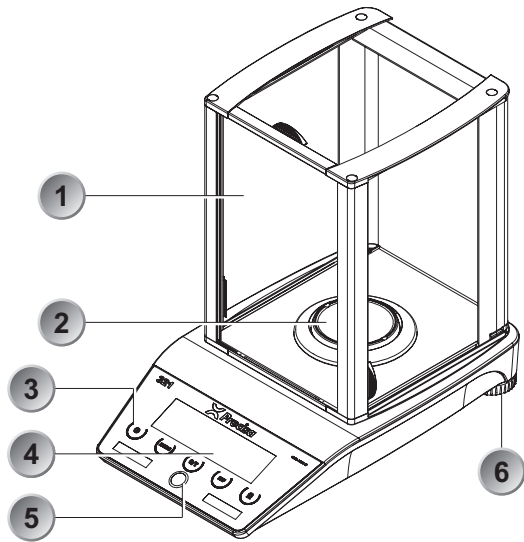
1.

Introduction

There are various models in the 321 Series. These brief instructions describe the balance with windshield. If you are using a less complex model, the descriptions of elements that are not included can be skipped. These brief instructions are intended as an initial overview and are an extract from the original operating instructions. The brief instructions must be available at all times at the workstation. More detailed descriptions can be found in the original operating instructions, which can be downloaded from our website. You should read the original operating instructions in full to ensure that you make the best possible use of your balance.

2.

Safety precautions

- When the balance is used in environments with increased safety requirements, the relevant precautions must be taken.
- The balance must only be operated with a fault-free power supply unit.
- Before connecting the power supply unit, ensure that the operating voltage specified on the power supply unit matches the mains supply.
- The balance must only be operated with the power supply unit supplied.
- If safe operation of the balance is not possible, it must be disconnected from the power supply and secured against unintentional use.

3.**Technical data**

| | |
|--|---|
| Casing dimensions W x D x H for: - Readability 0.1 mg - Readability 1 mg - Readability > 1 mg | 228 x 381 x 345 mm 228 x 381 x 124 mm 228 x 381 x 90 mm |
| Net weight: - Readability 0.1 mg - Readability 1 mg - Readability > 1 mg | 6.6 kg 5.5 kg 5.6 kg |
| Weighing range for each model (see details on nameplate) | 120 g to 10 200 g |
| Allowable ambient conditions (for calibratable balances, see details on the nameplate) | 0 °C to 40 °C |
| Allowable relative humidity | 25 % - 85 %, non-condensing |
| Mains connection (19) | 90 V - 264 V, 47 Hz - 63 Hz |
| Power consumption | 18 W, 12 V DC, 1.5 A |
| Data transfer connection (17 and 18) | RS 232/V24 interface, USB type B (device) |

Tab. 2
Technical data

4. Starting up

4.1 Choice of a suitable location

- Solid, firm and low-vibration, horizontal base
- No direct sunlight
- No draft
- Ambient temperature: 0 °C - 40 °C
- Relative humidity (non-condensing): 25 % - 85 %

4.2 Unpacking the balance

NOTE

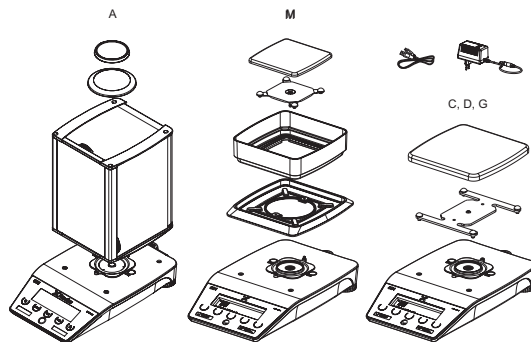
Retain the original packaging in order to avoid damaging the balance during shipping or transport and to enable the balance to be stored in ideal conditions when out of use for long periods.

- Carefully unpack the balance.

4.3

Assembling the balance**CAUTION**

Before connecting the device to the mains, be sure that the operating voltage stated on the device and on the power adaptor agrees with the local mains voltage. Do not touch any key during the following start-up sequence until after about 3 seconds an acoustic signal sounds.


- Place the dust protection cover on the balance.
- Install the draft shield (1).
- Place the balance pan holder in position and add the weighing pan (2).
- Insert the supplied power supply unit into the socket (19) on the back of the balance.

4.4 Under-floor weighing



For weighing objects which, because of their size or shape, cannot be put in or on the balance.

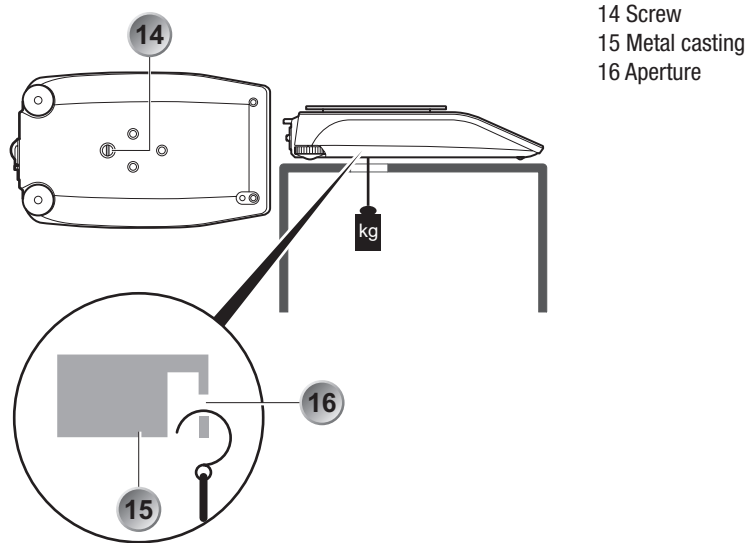- Remove the screw (14) on the underside of the balance.
- Place the balance over an opening.
- Hang the hook for under-bench weighing, to which the object can be secured, into the aperture (16) of the metal casting (15).


4.5 Levelling the balance

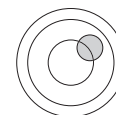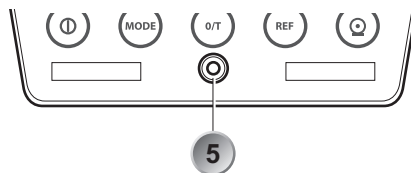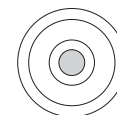- Adjust the two adjustable feet (6) so that the air bubble in the level control (5) below the buttons on the control panel (3) lies precisely in the centre of the sight glass marking.

Wrong

Right

4.6 Calibrating the balance

- Press the  key to switch on the balance.
- Press the  key until the 20-character info line (12) in the display (4) shows «CALIBRATION».

The balance begins the calibration process.

- Follow the instructions in the info line (12) and the eight-digit measurement display (13).

NOTE

The balance must be recalibrated each time its location is changed.

5. Key functions

| Key | Designation | Function | Position index |
|---|-------------|---|----------------|
|  | ON/OFF | Switches the balance on/off | 7 |
|  | MODE | Menu or change to an application / change to normal weighing | 8 |
|  | O/T | Activates the tare functions and calibration function and other functions | 9 |
|  | REF | Activates a function that corresponds to the application currently in use | 10 |
|  | PRINT | Activates printing functions and the Back button in the menu | 11 |

Tab. 5
Key functions

6.

Overview of menus

**NOTE**

With user guidance activated, the keys that currently have a function or can be used backlit.

6.1

Menu control

Navigating through menu levels:

| Key | Press duration | Function |
|---|----------------|---|
|  | short | Switch to a sub-menu or start an input |
|  | long | Switch to another menu item in the same menu |
|  | short | Switch from the sub-menu to the main menu or exit the main menu |

Tab. 6.1.1
Navigating through
menu levels

Make an input:

| Key | Press duration | Function |
|---|----------------|---|
|  | short | Change the value or digit |
|  | long | Change the position or end the input or |
|  | short | End the input |

Tab. 6.1.2
Making an input

6.2

Configuration menu

- Switch the balance off.
- First press the  key briefly and then immediately press the  key with a long press:

| Menu items | Function |
|-------------------|---|
| LANGUAGE | Selecting the language |
| UNIT-1 | Unit in which the weighing results are displayed |
| SET DATA PRINT | Print formats, type of values to be printed out |
| SET CALIBRATION | Calibration method |
| SET WEIGHING MODE | Stability mode, Auto-Standby mode, zero correction, tare method, zero range |
| SET INTERFACE | Baud rate, parity, handshake functions of the peripheral interface |
| SET DATE AND TIME | Date and time |
| PASSWORD | Password protection for menu definitions |
| THEFTCODE | Activating / deactivating and changing the anti-theft code |
| KEY TONE | Keypad sound |
| BUS | Enable or disable the Precisa BUS |

Tab. 6.2
Contents of the
configuration menu

6.3

Application menu

- Check that the balance is switched on.
- Press the  key with a long press:

| Menu items | Function |
|-------------------|--|
| SET APP. UNITS | Select an application |
| SETUP APPLICATION | Specify the parameters for the application selected under "APPLICATION" |
| AUTO-START ON | The selected application can, if required, be loaded automatically every time the balance is switched on |

Tab. 6.3
Contents of the
application menu

7. Key menus

7.1 Tare menu

7.1.1 Taring the balance

- Press the  key with a short press:
→ The balance is tared.

7.1.2 Other functions

- Press the  key until you reach the required function:

| Menu items | Function |
|-------------|---|
| CALIBRATION | The balance begins the calibration process. |

- Release the  key.

Tab. 7.1.2
Contents of the
tare menu

7.2 Mode menu

7.2.1 Switch to an activated application or back to weighing mode

- Press the  key with a short press:
→ The balance switches to the last application activated. This acts like a switchover.

7.2.2 Other functions

- Press the  key until you reach the required function:

| Menu items | Function |
|---------------------|--|
| WEIGHING | Switches to weighing mode |
| UNITS (COUNT; etc.) | Switches to the selected application units, counting, etc. |
| AUTO-START ON | The selected application can, if required, be loaded automatically every time the balance is switched on |

- Release the  key.

Tab. 7.2.2
Contents of the
mode menu

7.3

Print menu

7.3.1

Print individual value / report

- Press the  key with a short press:
→ The individual value or report is printed out.

7.3.2

Other functions

- Press the  key until you reach the required function:

| Menu items | Function |
|---------------------|---------------------------------------|
| RESET PROD. COUNTER | Reset the product counter to 1 |
| PRINT STATUS | Print out balance settings |
| PRINT APPLICATIONS | Print out application settings |
| PRINT CALIBRATIONS | Print out calibration information |
| PRINT FIRMWARE HIST | Print out firmware update information |

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