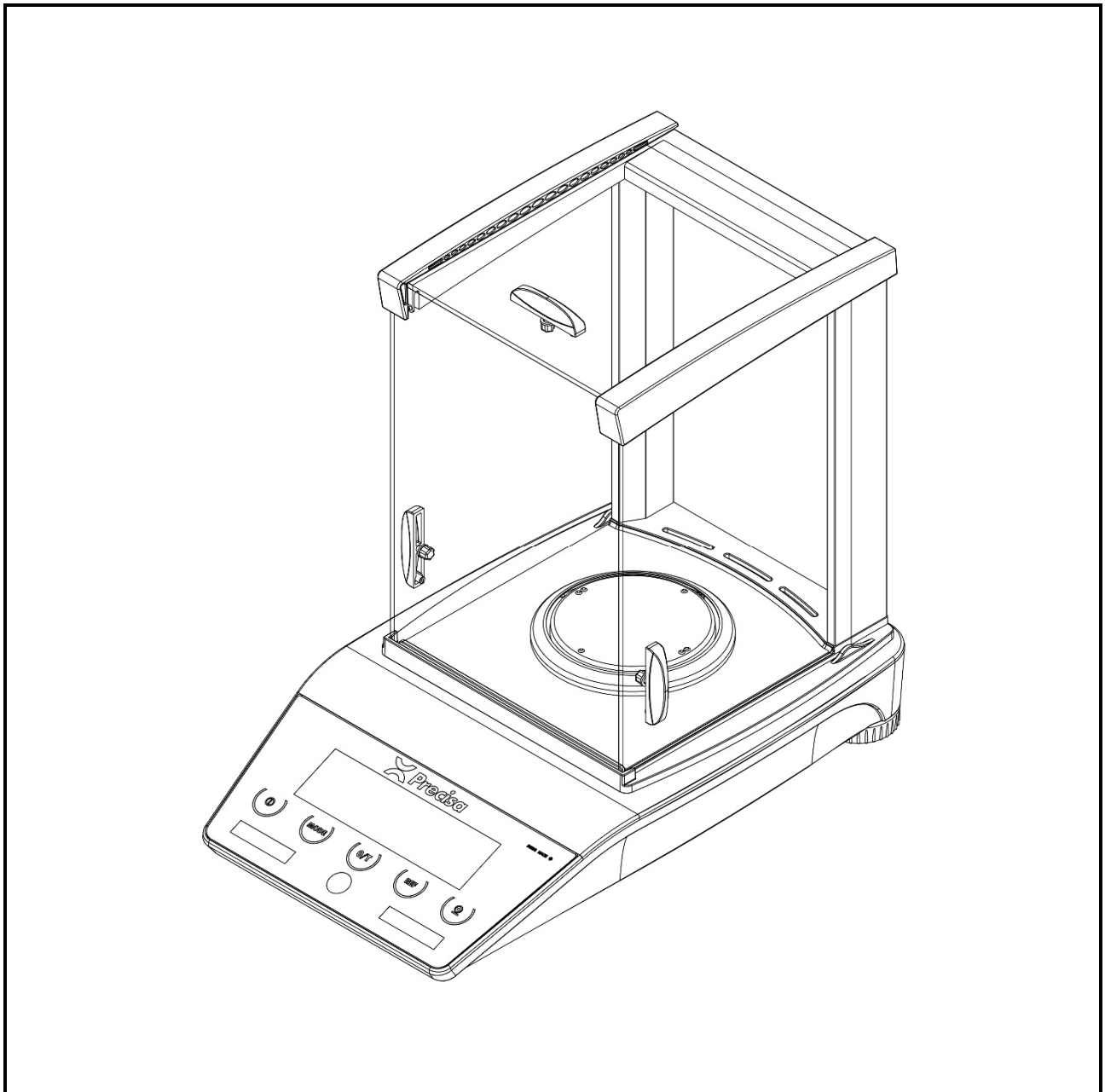
SWISS MADE

# 321 LS

Quick Start  
Schnelleinstieg  
Départ rapide  
Inicio rápido

M350-8173-900d1

 **Precisa**  
THE BALANCE OF QUALITY

## Safety precautions

When the balance is used in environments with increased safety requirements, the relevant precautions must be taken.

The balance must only be operated with a fault-free power adapter.

Before connecting the power adapter, ensure that the operating voltage specified on the power adapter matches the mains supply.

The balance must only be operated with the power adapter supplied.

If safe operation of the balance is not possible, it must be disconnected from the power supply and secured against unintentional use.

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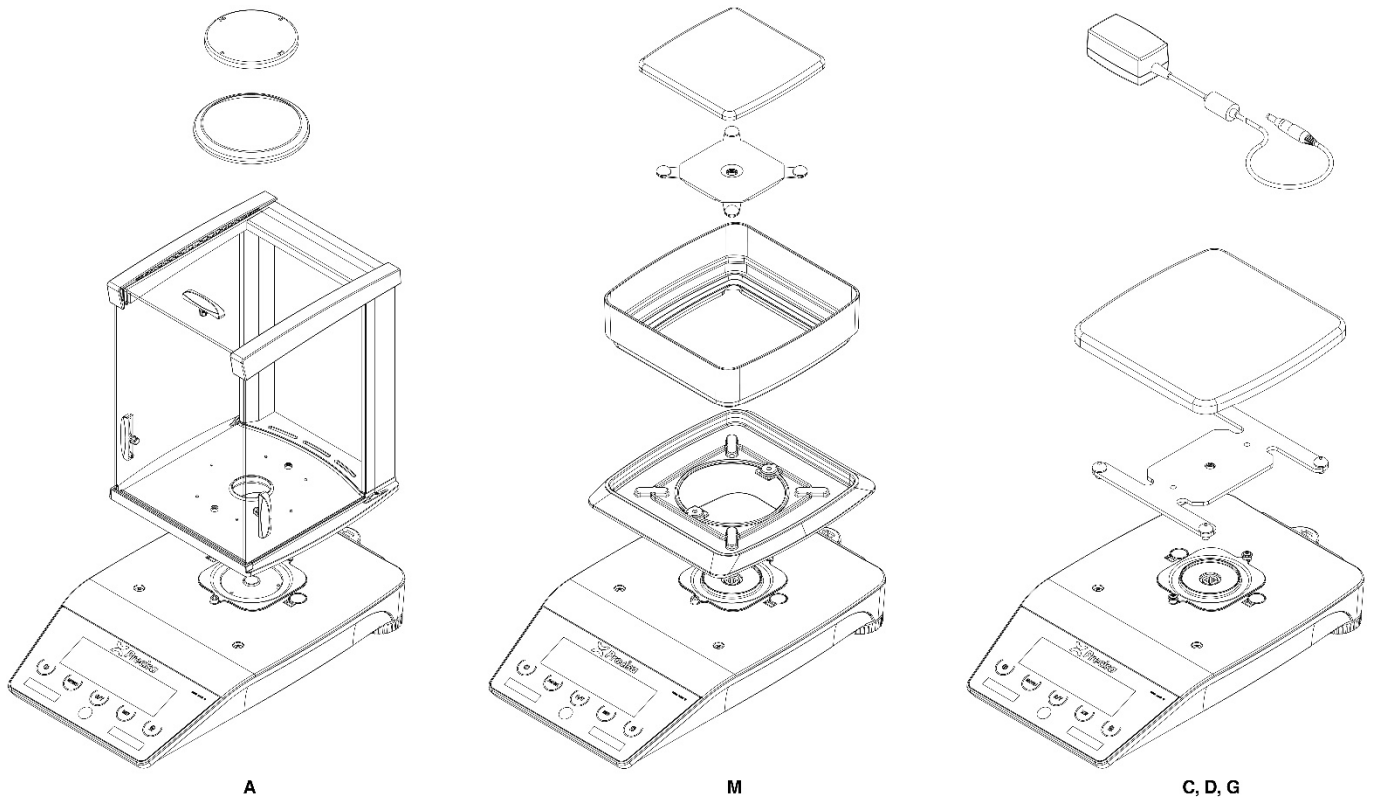
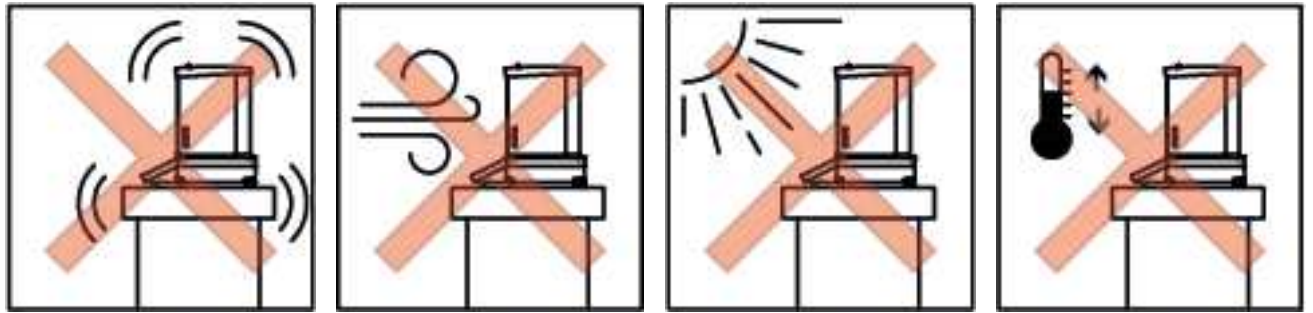
## Starting up – Inbetriebnahme – Mise en service – Puesta en servicio


Only work with a balance at operating temperature (warm-up time at least 2 hours). If you have not used the balance for a long time, briefly load it with your hand and tare it.

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## Operation

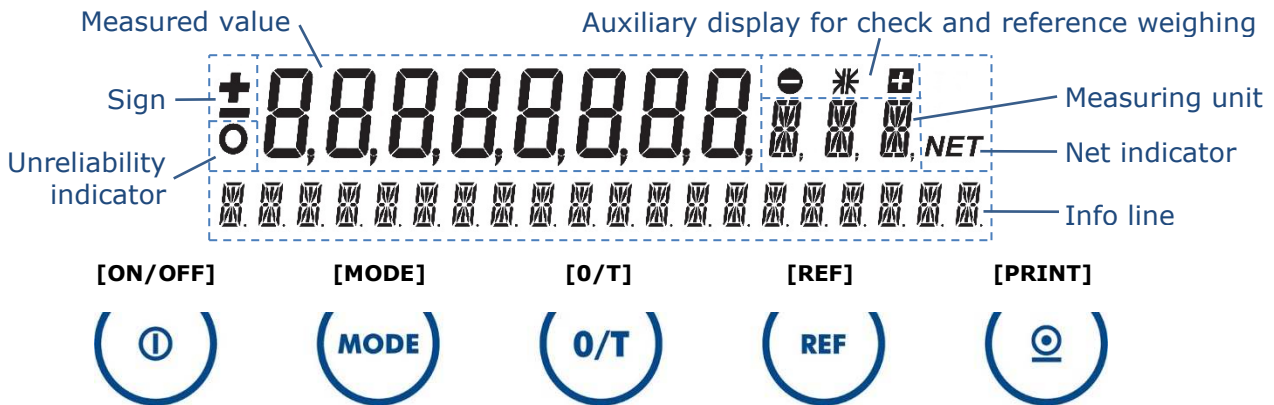The measured value is stable as soon as the unit, e.g. "g", appears. The unreliability indicator lights up while the scale is recalculating the reference weight, for example, during counting, or when it is verified and in the warm-up phase.

### Application example: Adding up components (application net total)

If the balance is not set to English: Enter the configuration menu, switch the balance off and on with **[ON/OFF]** and, as soon as all segments of the display light up, press and hold **[MODE]** until "SPRACHE DEUTSCH/LANGUE FRANÇAISE" appears. Press **[MODE]**; "SPRACHE DEUTSCH/LANGUE FRANÇAISE" starts flashing. Press **[MODE]** repeatedly until "LANGUAGE ENGLISH" appears and confirm with **[PRINT]**; "LANGUAGE ENGLISH" stops flashing. Press **[PRINT]**; the balance finishes starting up.

To enter the application menu, keep **[MODE]** pressed until "SET APP: OFF" appears. Press **[MODE]**; "OFF" starts flashing. Press **[Down]** repeatedly until "NET-TOTAL" appears and confirm with **[PRINT]**. Keep **[MODE]** pressed until "SETUP APPLICATION" appears. Keep **[MODE]** pressed until "AUTO-START ON" appears. For manual value transfer press **[MODE]**; "ON" starts flashing, press **[MODE]**; "OFF" appears, confirm with **[PRINT]**; "OFF" stops flashing. Press **[PRINT]**; the balance exits the application menu. Press **[MODE]**; the balance starts the application Net total and displays 0g as the measured value and "STO 0 RES INF" in the info line.

Place the first component, e.g. 50g, and press **[REF]**; the balance records the value, tares itself and displays the capacity of the weighing range analogically in the info line. Place the second component, e.g. 100g, and press **[REF]**, and finally place the third component, e.g. 20g, and press **[REF]**.

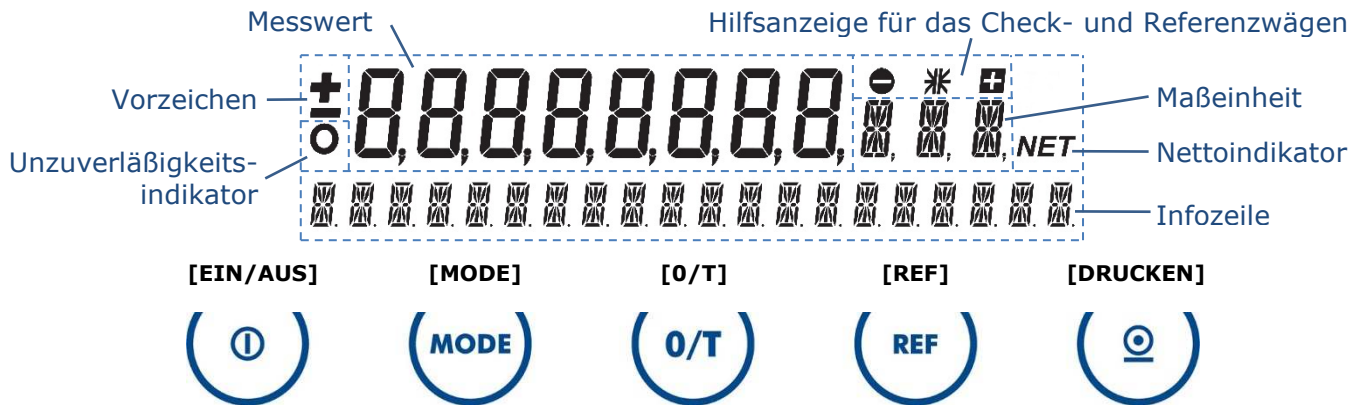
While adding up, the current total weight can be viewed at any time by pressing **[MODE]**. The balance then switches to weighing mode. To return to the application, press **[MODE]** again. To switch from the analogue display to the application commands in the info line, press **[MODE]** twice; the balance switches to the weighing mode and back to the application. It then displays e.g. "STO 3 RES INF".

Press and hold **[REF]** until "INF" starts flashing; the balance displays in the info line all information on the weights just added up. Use **[REF]** to scroll through the entire stack of info lines: Total, remaining capacity and all individual weights, e.g. 50g, 100g and 20g. Press **[MODE]** to exit the information query; the balance switches back to the running application with analogue display and is ready for another weighing-in.

To restart, press and hold **[REF]** until "RES" starts flashing; the balance remains in the application with analogue display and resets the counter to 0.

To quit the application, press **[MODE]**; the balance switches back to weighing mode.

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© Precisa Gravimetrics AG, 8953 Dietikon, Switzerland.