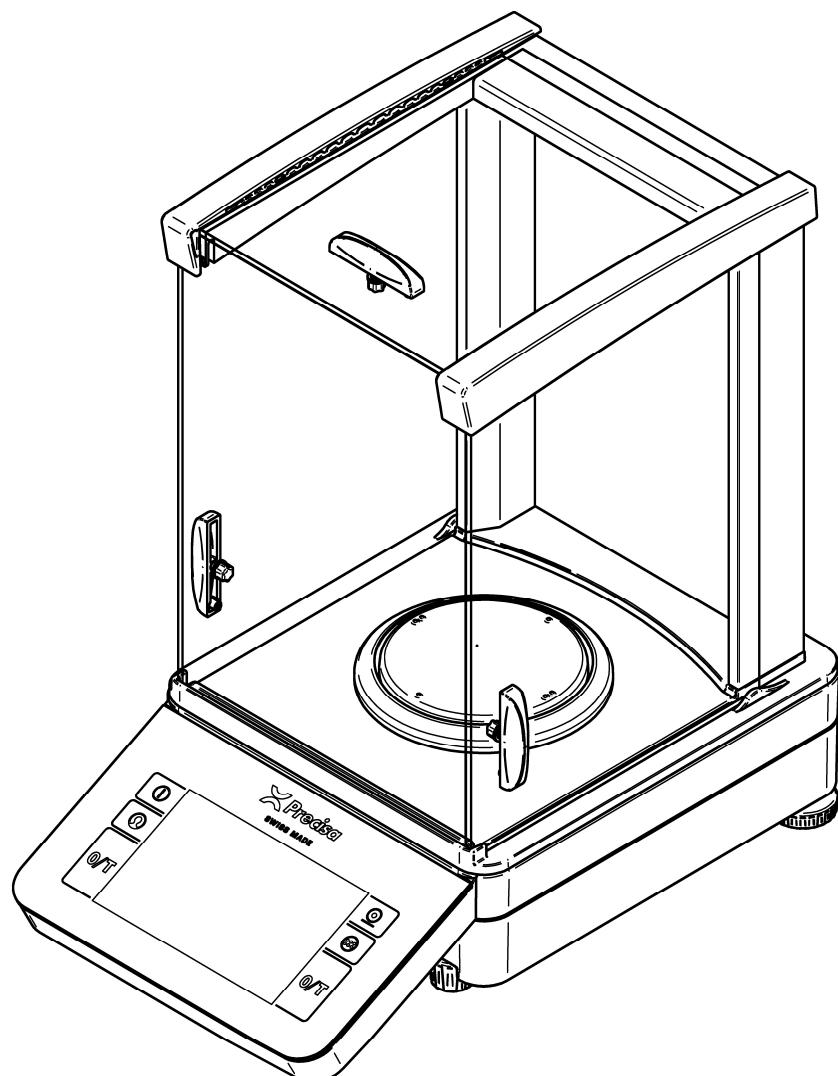
SWISS MADE

# 520 PT

Quick Start  
Schnelleinstieg  
Départ rapide  
Inicio rápido

M350-8186-900a1

 **Precisa**  
THE BALANCE OF QUALITY

## Safety precautions

When the balance is used in environments with increased safety requirements, the relevant precautions must be taken.

The balance must only be operated with a fault-free power adapter.

Before connecting the power adapter, ensure that the operating voltage specified on the power adapter matches the mains supply.

The balance must only be operated with the power adapter supplied.

If safe operation of the balance is not possible, it must be disconnected from the power supply and secured against unintentional use.

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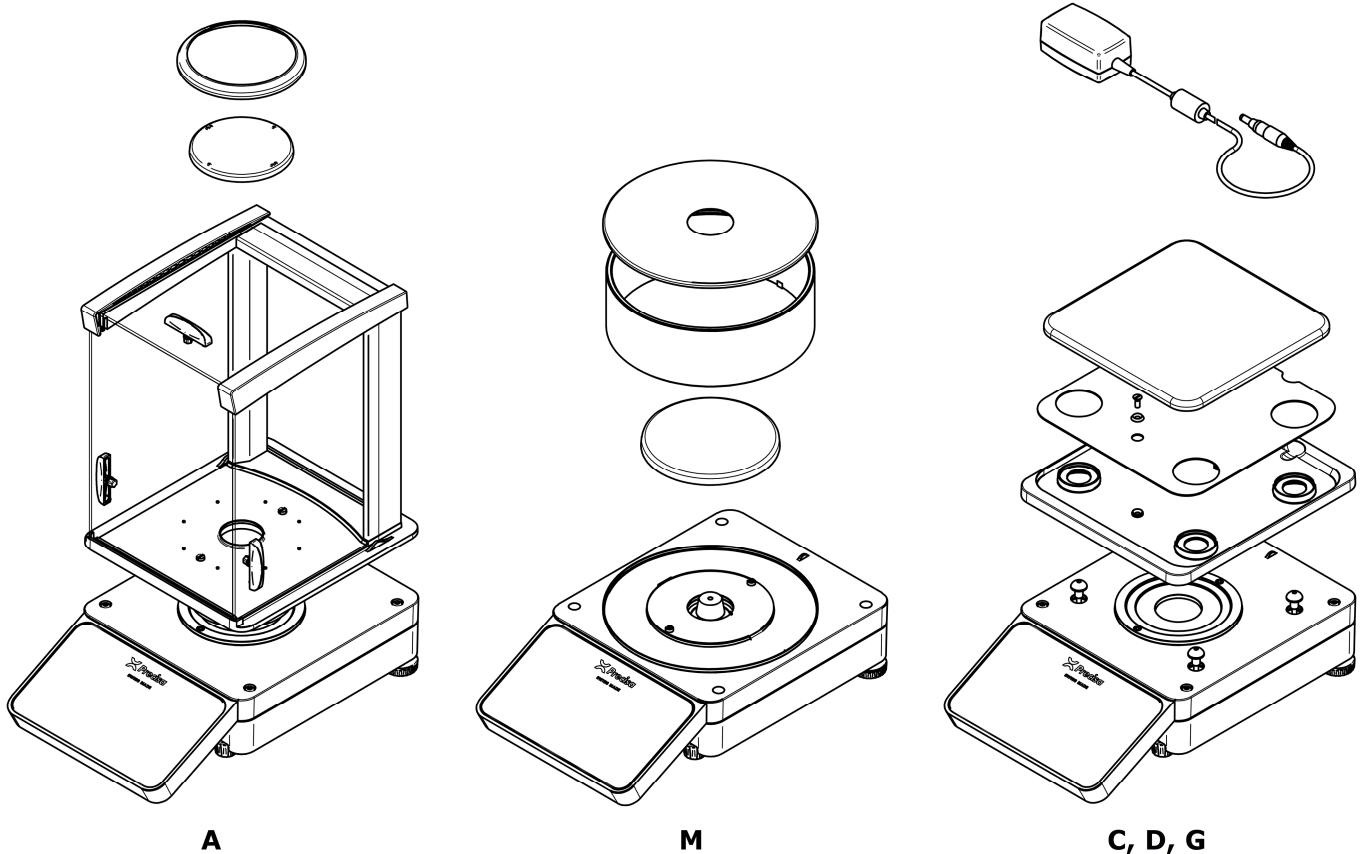
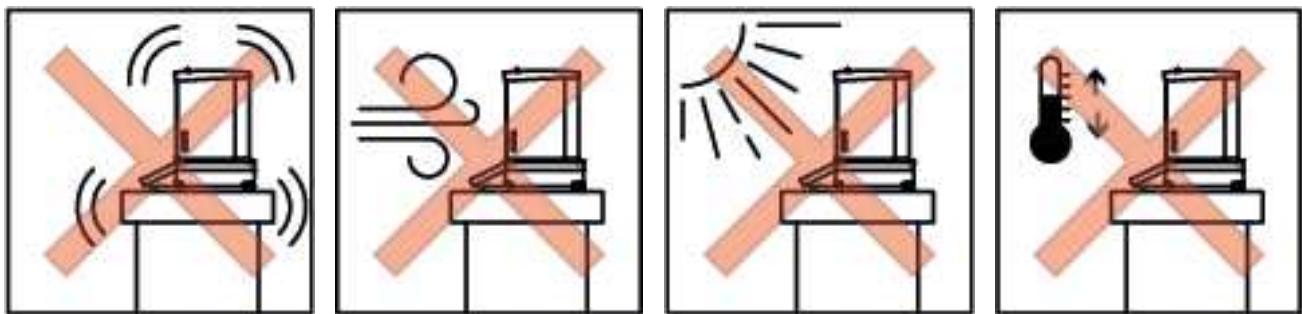
## Starting up – Inbetriebnahme – Mise en service – Puesta en servicio










Only work with a balance at operating temperature (warm-up time at least 2 hours). If you have not used the balance for a long time, briefly load it with your hand and tare it.

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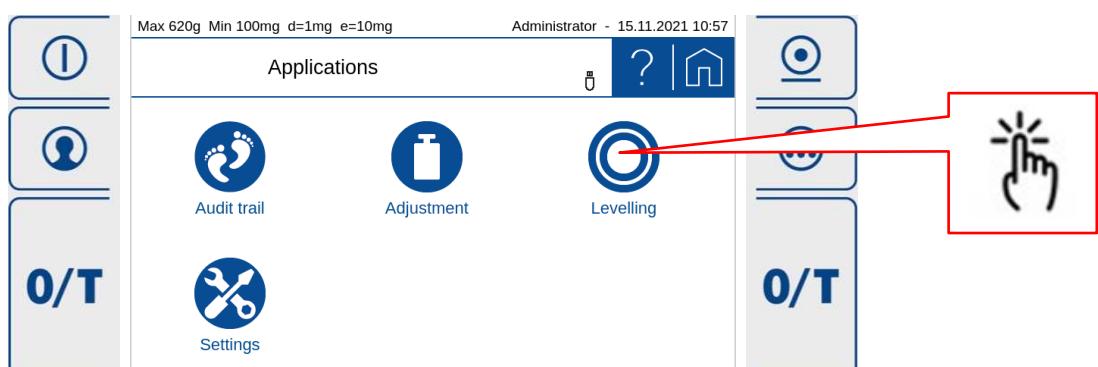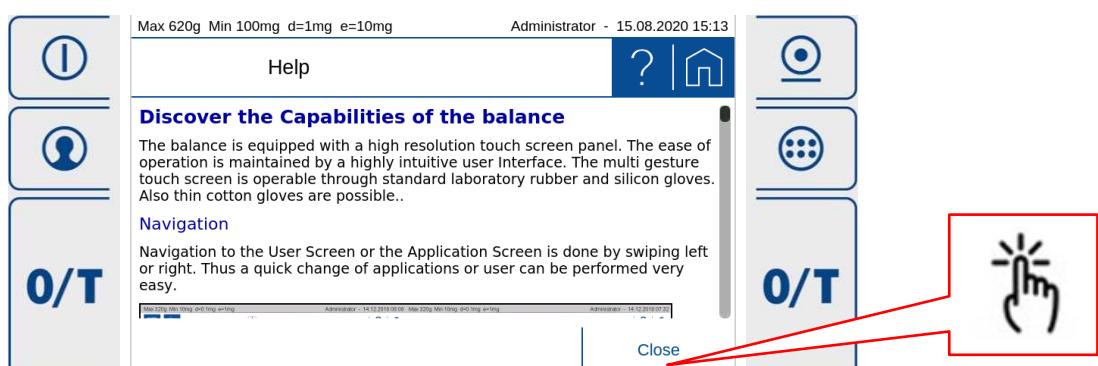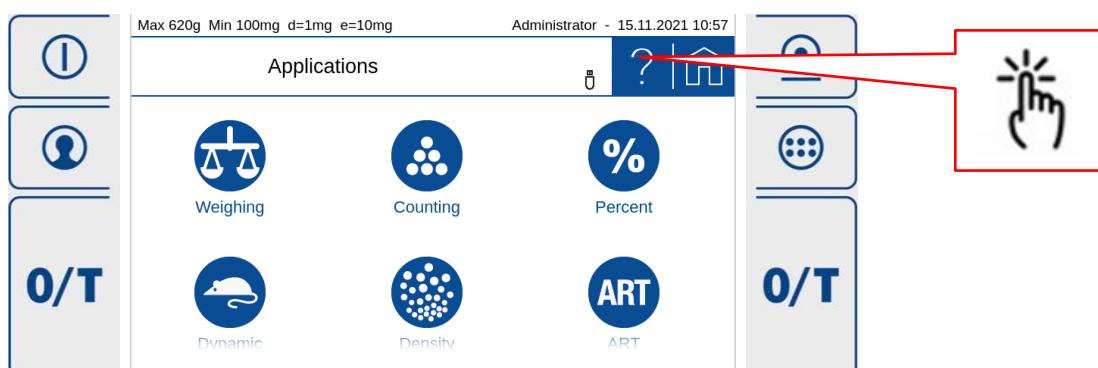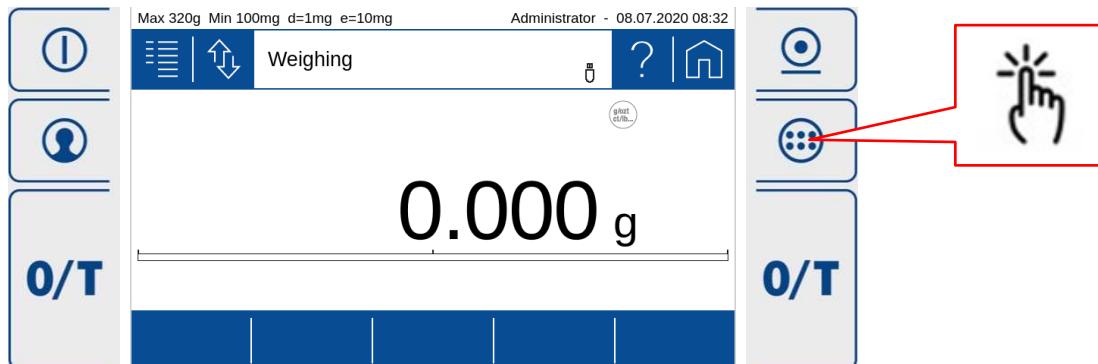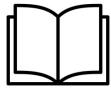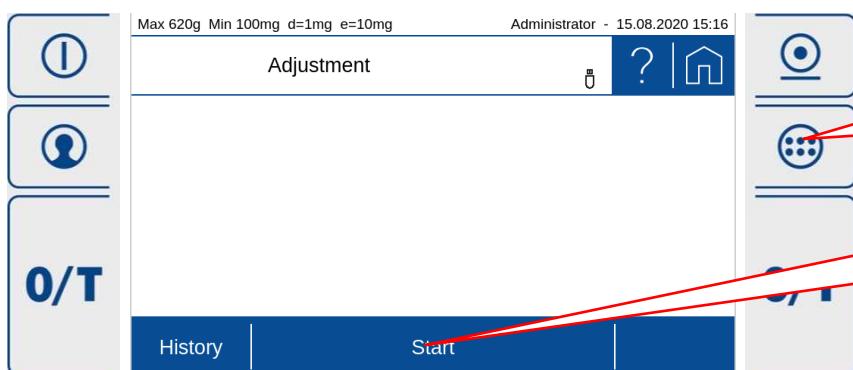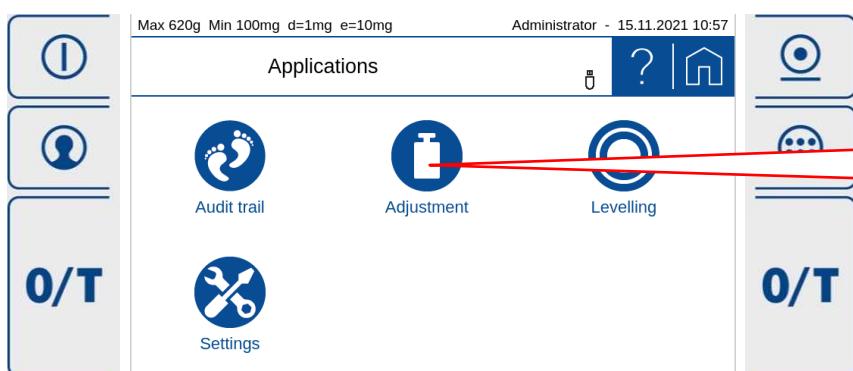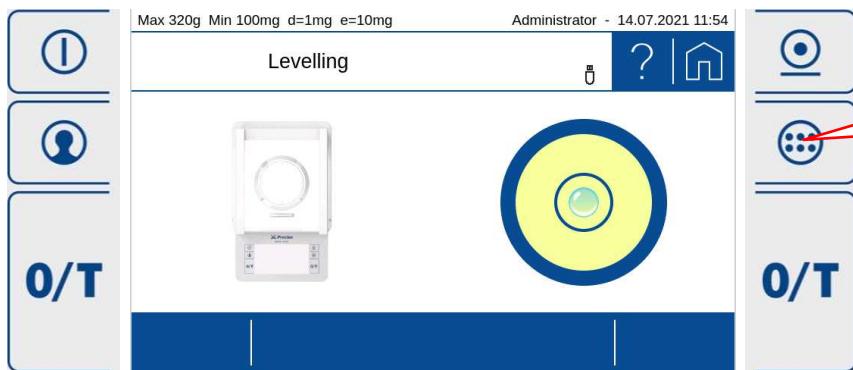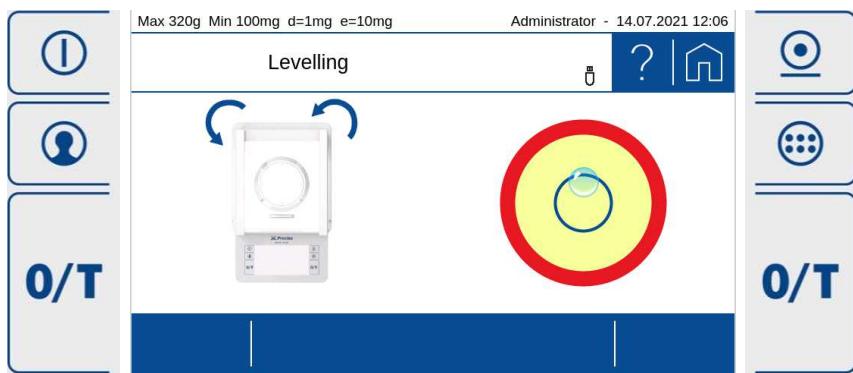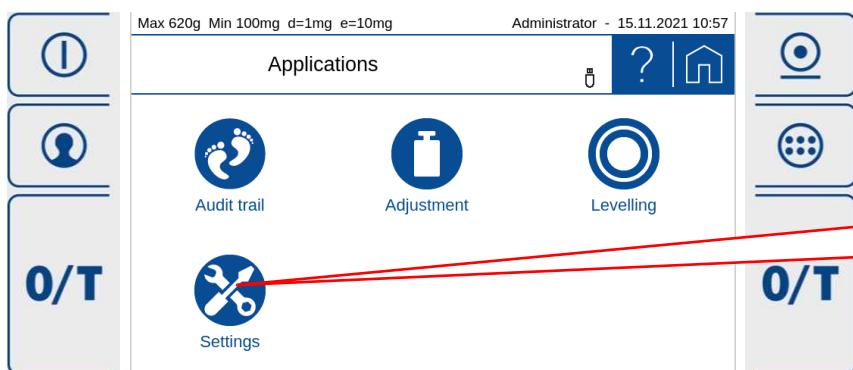
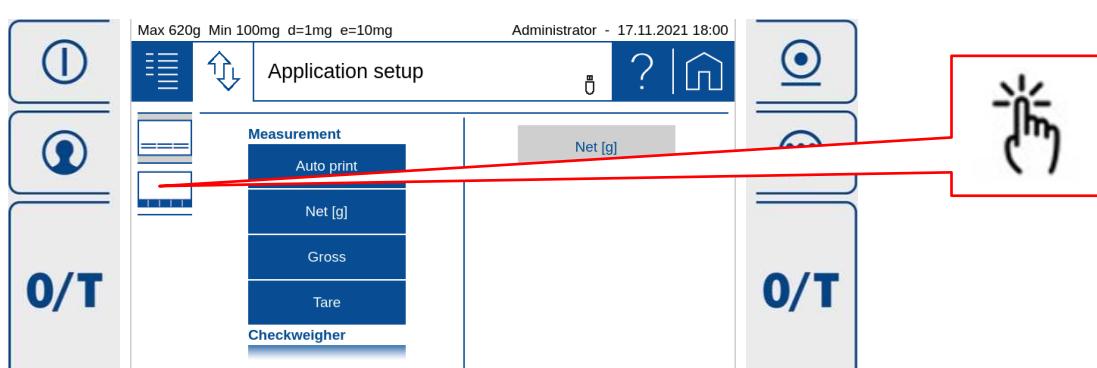
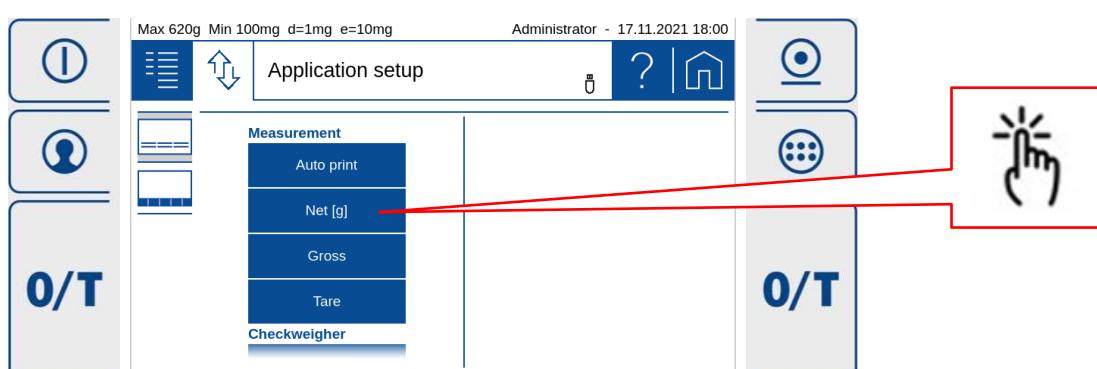
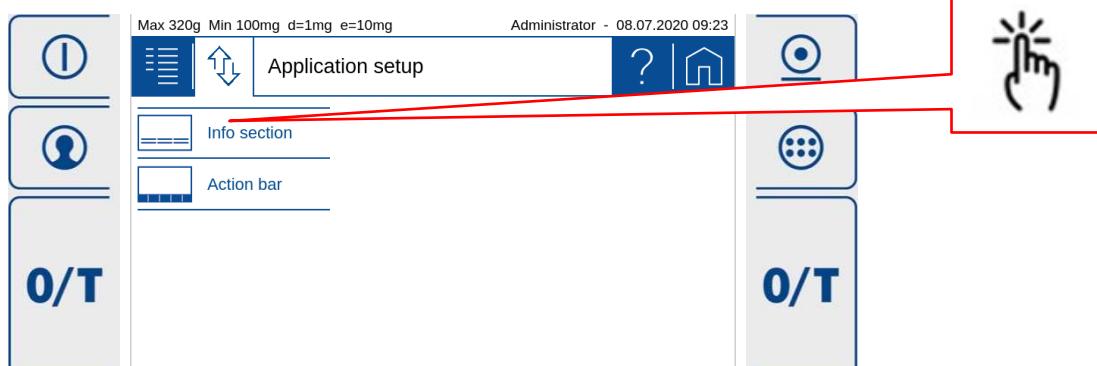
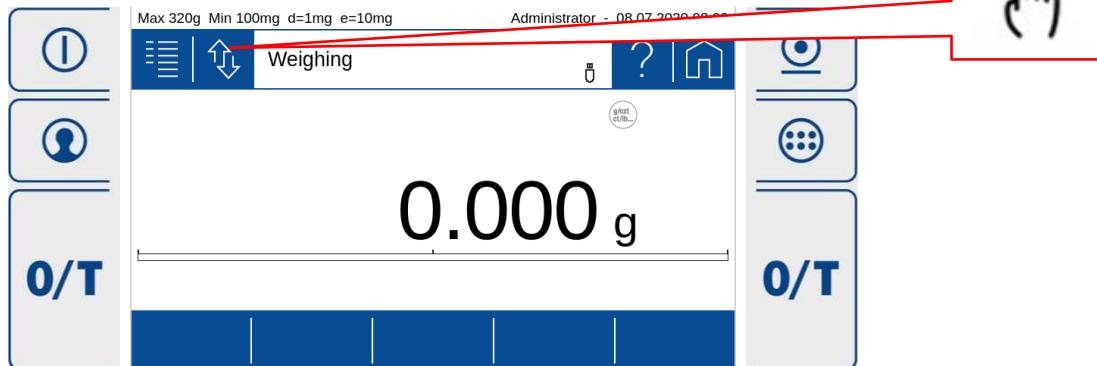
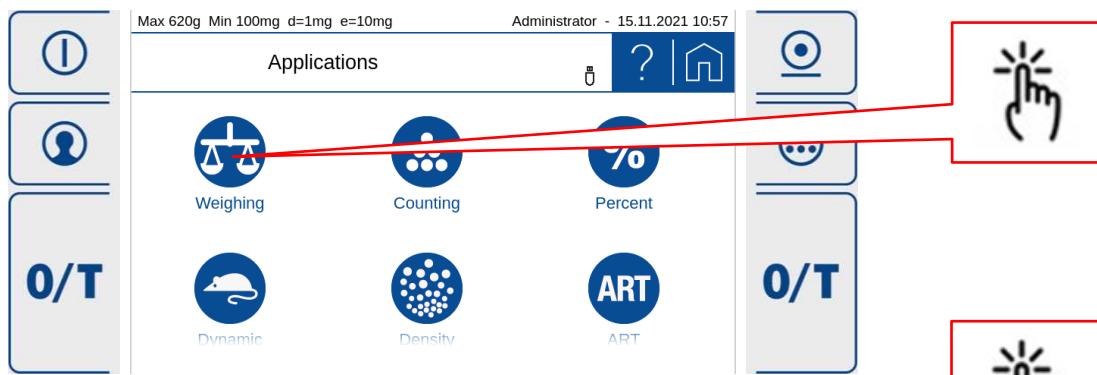
1.

**Top Left Panel:**

- Icons:** Information (I), User (User icon), Off/On (O/T).
- Text:** Max 620g Min 100mg d=1mg e=10mg, Administrator - 17.11.2021 18:03.
- Buttons:** Application setup, Store 1, Show statistics, Clear statistics, Delete last sample, Statistics, Print stat, Print chart.
- Display:** mg
- Right Side:** A red box highlights the 'Application setup' button, which is connected by a red line to a hand icon pointing at it.
  

**Second Row Left Panel:**

  - Icons:** Information (I), User (User icon), Off/On (O/T).
  - Text:** Max 620g Min 100mg d=1mg e=10mg, Administrator - 17.11.2021 18:05.
  - Buttons:** Application setup, Store 1, Show statistics, Clear statistics, Delete last sample, Statistics, Print stat, Print chart.
  - Display:** mg
  - Right Side:** A red box highlights the 'Application setup' button, which is connected by a red line to a hand icon pointing at it.
  

**Third Row Left Panel:**

    - Icons:** Information (I), User (User icon), Off/On (O/T).
    - Text:** Max 620g Min 100mg d=1mg e=10mg, Administrator - 17.11.2021 18:05.
    - Buttons:** Application setup, Store 1, Show statistics, Clear statistics, Delete last sample, Statistics, Print stat, Print chart.
    - Display:** mg
    - Right Side:** A red box highlights the 'Application setup' button, which is connected by a red line to a hand icon pointing at it.
  

**Fourth Row Left Panel:**

      - Icons:** Information (I), User (User icon), Off/On (O/T).
      - Text:** Max 620g Min 100mg d=1mg e=10mg, Administrator - 19.08.2020 14:00.
      - Buttons:** Application setup, Store 1, Show statistics, Clear statistics, Delete last sample, Statistics, Print stat, Print chart.
      - Display:** mg
      - Right Side:** A red box highlights the 'Application setup' button, which is connected by a red line to a hand icon pointing at it.
  

**Fifth Row Left Panel:**

        - Icons:** Information (I), User (User icon), Off/On (O/T).
        - Text:** Max 620g Min 100mg d=1mg e=10mg, Administrator - 19.08.2020 14:01.
        - Buttons:** Application setup, Store 1, Show statistics, Clear statistics, Delete last sample, Statistics, Print stat, Print chart.
        - Display:** Net [g] 0.000 g
        - Right Side:** A red box highlights the 'Net [g]' field, which is connected by a red line to a hand icon pointing at it.
  

**Sixth Row Left Panel:**

          - Icons:** Information (I), User (User icon), Off/On (O/T).
          - Text:** Max 620g Min 100mg d=1mg e=10mg, Administrator - 19.08.2020 14:01.
          - Buttons:** Application setup, Store 1, Show statistics, Clear statistics, Delete last sample, Statistics, Print stat, Print chart.
          - Display:** NET 2556 mg
          - Right Side:** A red box highlights the 'Net [g]' field, which is connected by a red line to a hand icon pointing at it.

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18 VA, 12 V DC, 1.5 A

## Note – Hinweis – Remarque – Nota

Retain the original packaging in order to avoid damaging the balance during shipping or transport and to enable the balance to be stored in ideal conditions (5 .. 40°C / 41 .. 104°F) when out of use for long periods.

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No reproduction in any form of this document, in whole or in part, may be made without written authorization from Precisa Gravimetrics AG.

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© Precisa Gravimetrics AG, 8953 Dietikon, Switzerland