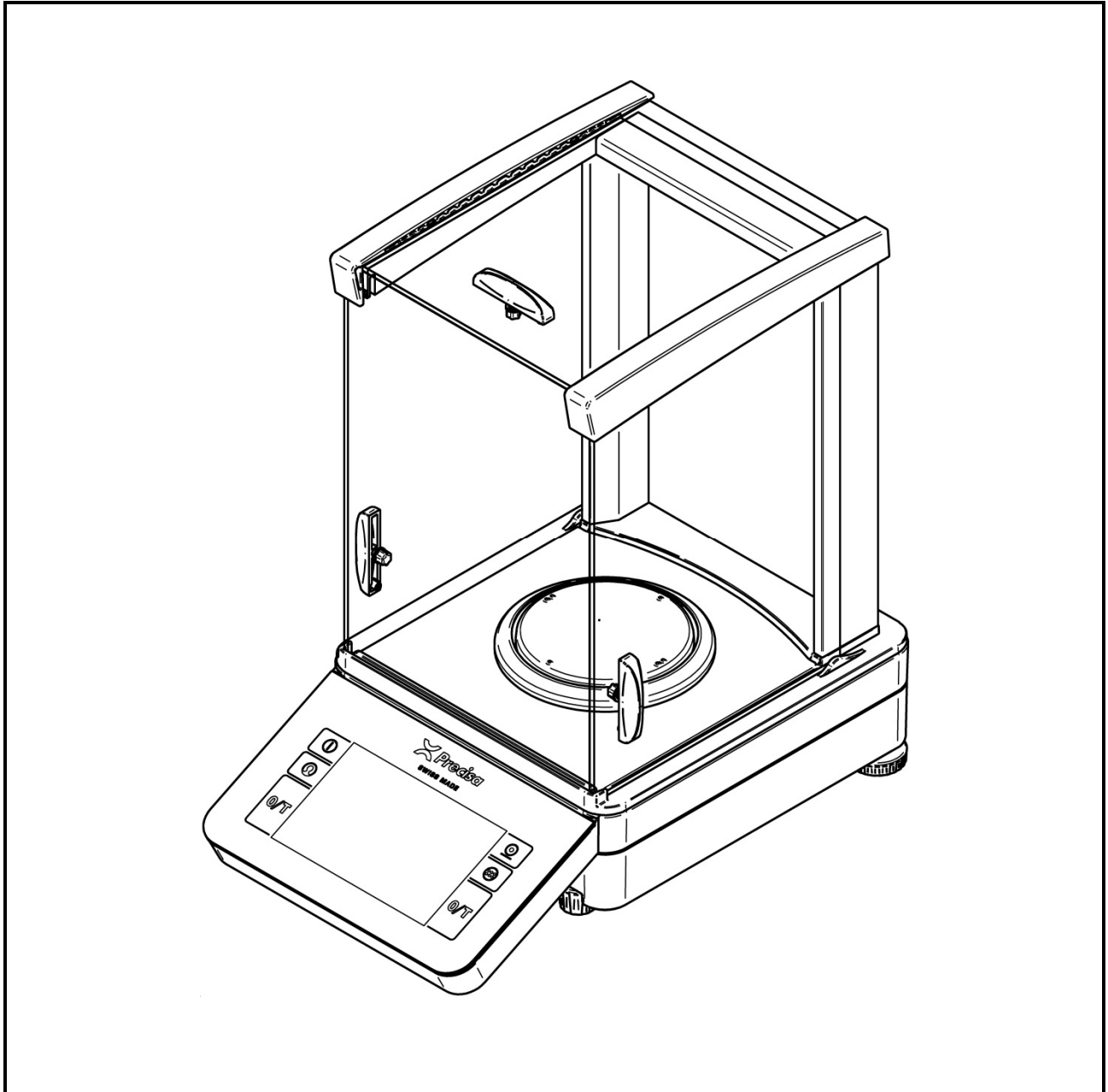
SWISS MADE

520 PT

Quick Start
Schnelleinstieg
Mode d'emploi simplifié
Inicio rápido

M350-8186-200a1

 **Precisa**

Safety precautions

When the balance is used in environments with increased safety requirements, the relevant precautions must be taken.

The balance must only be operated with a fault-free power adapter.

Before connecting the power adapter, ensure that the operating voltage specified on the power adapter matches the mains supply.

The balance must only be operated with the power adapter supplied.

If safe operation of the balance is not possible, it must be disconnected from the power supply and secured against unintentional use.

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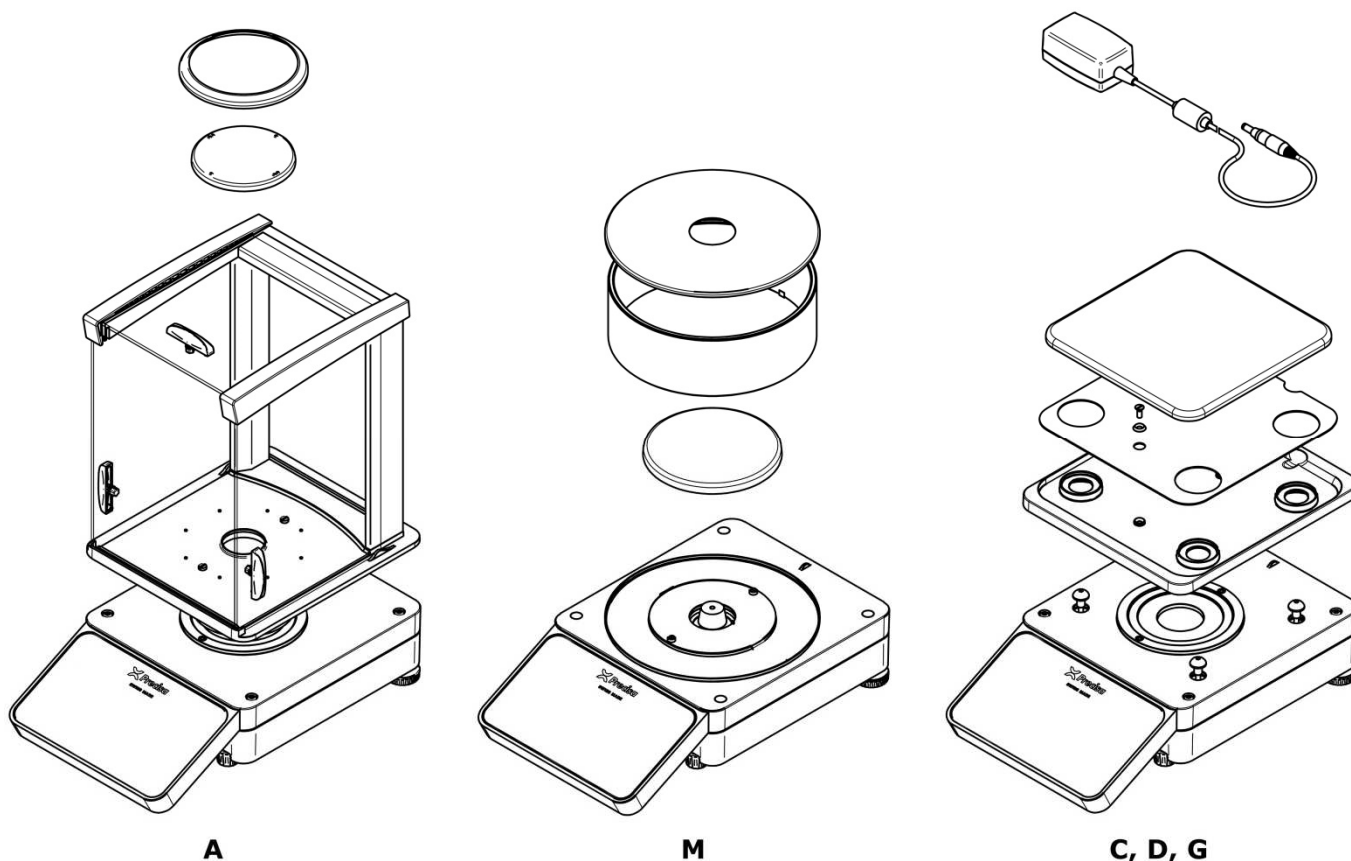
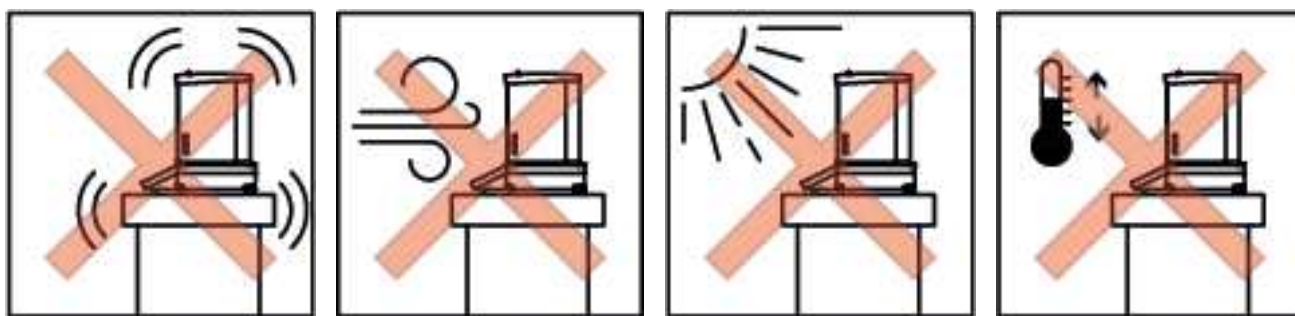
Starting up – Inbetriebnahme – Mise en service – Puesta en servicio

Only work with a balance at operating temperature (warm-up time at least 2 hours). If you have not used the balance for a long time, briefly load it with your hand and tare it.

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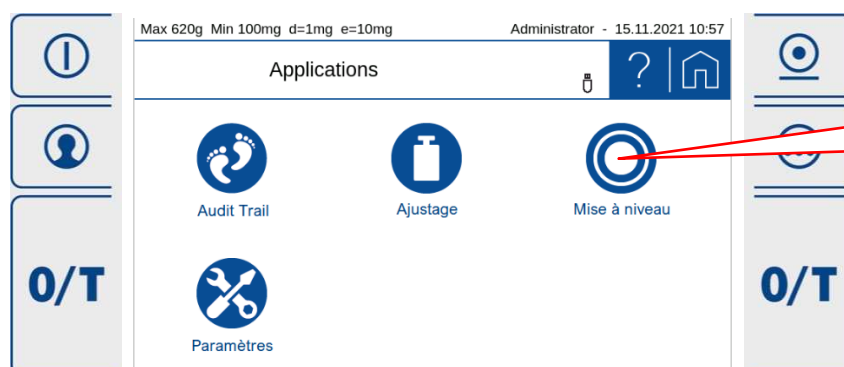
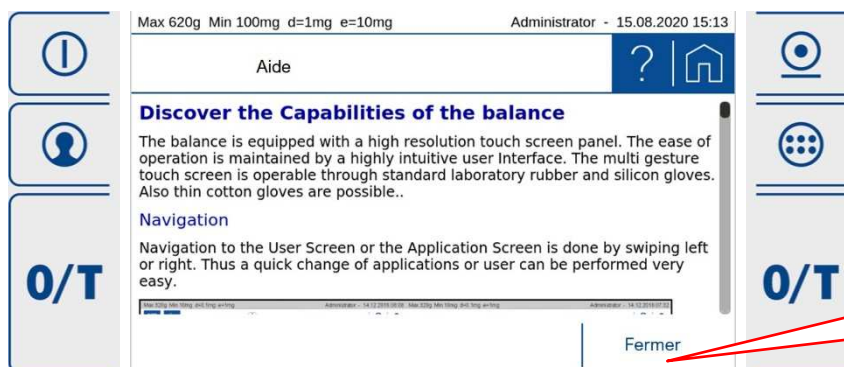
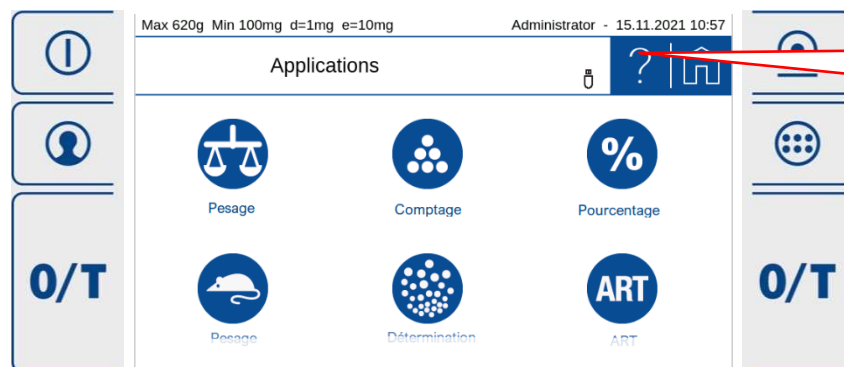
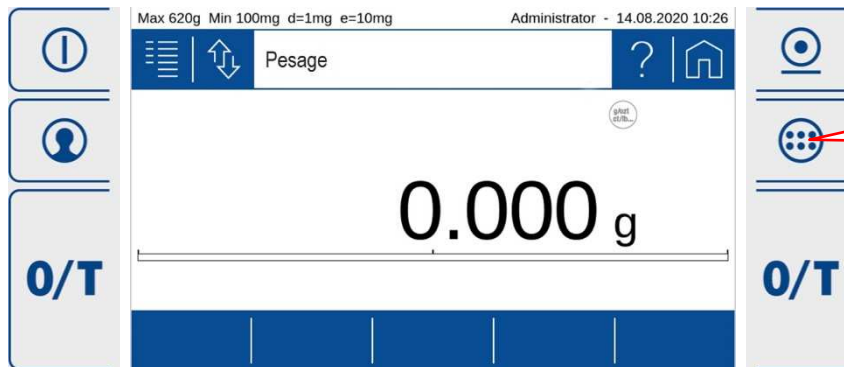
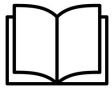
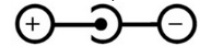
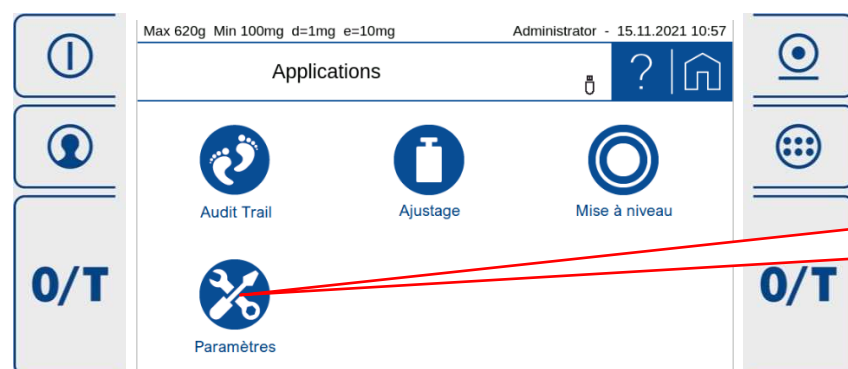
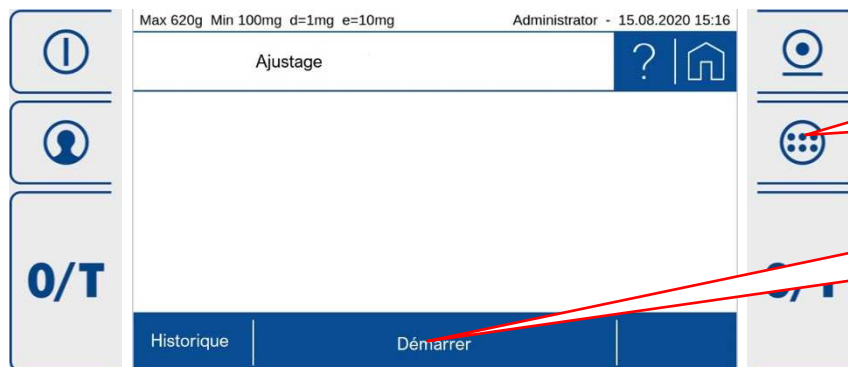
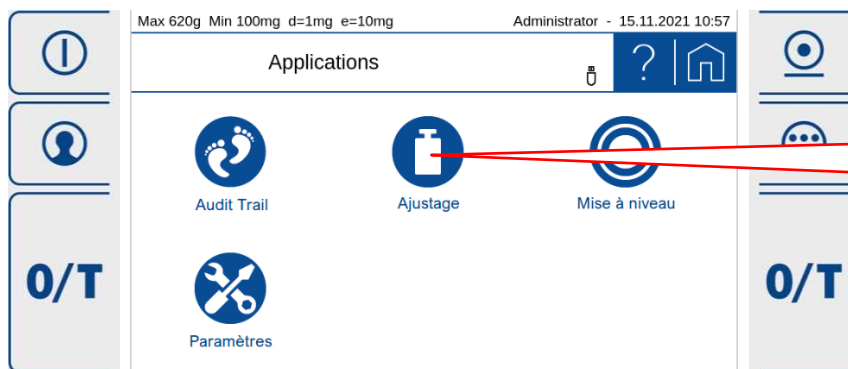
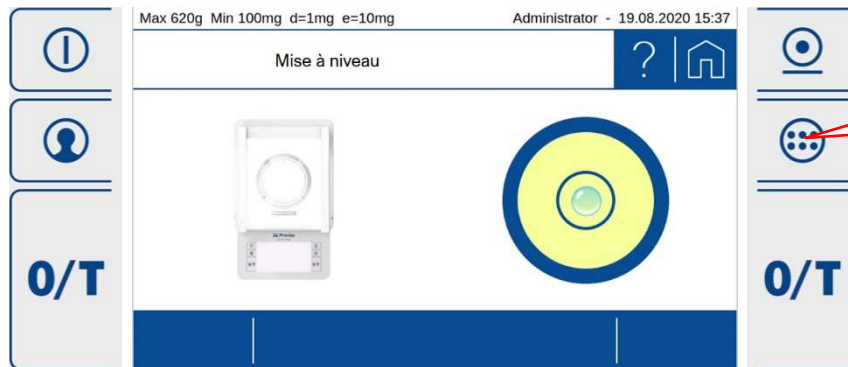
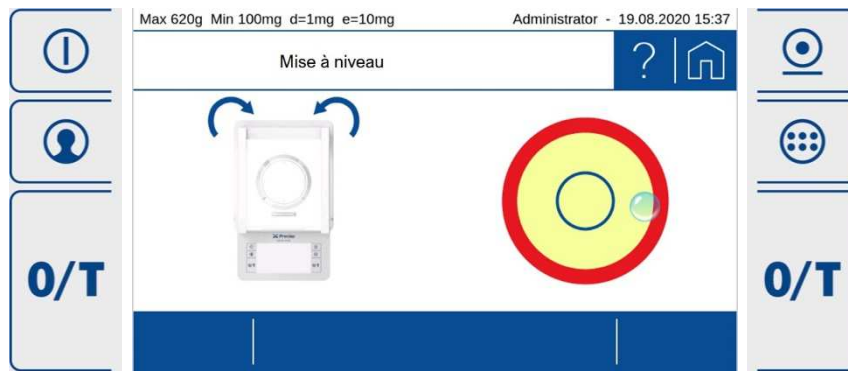
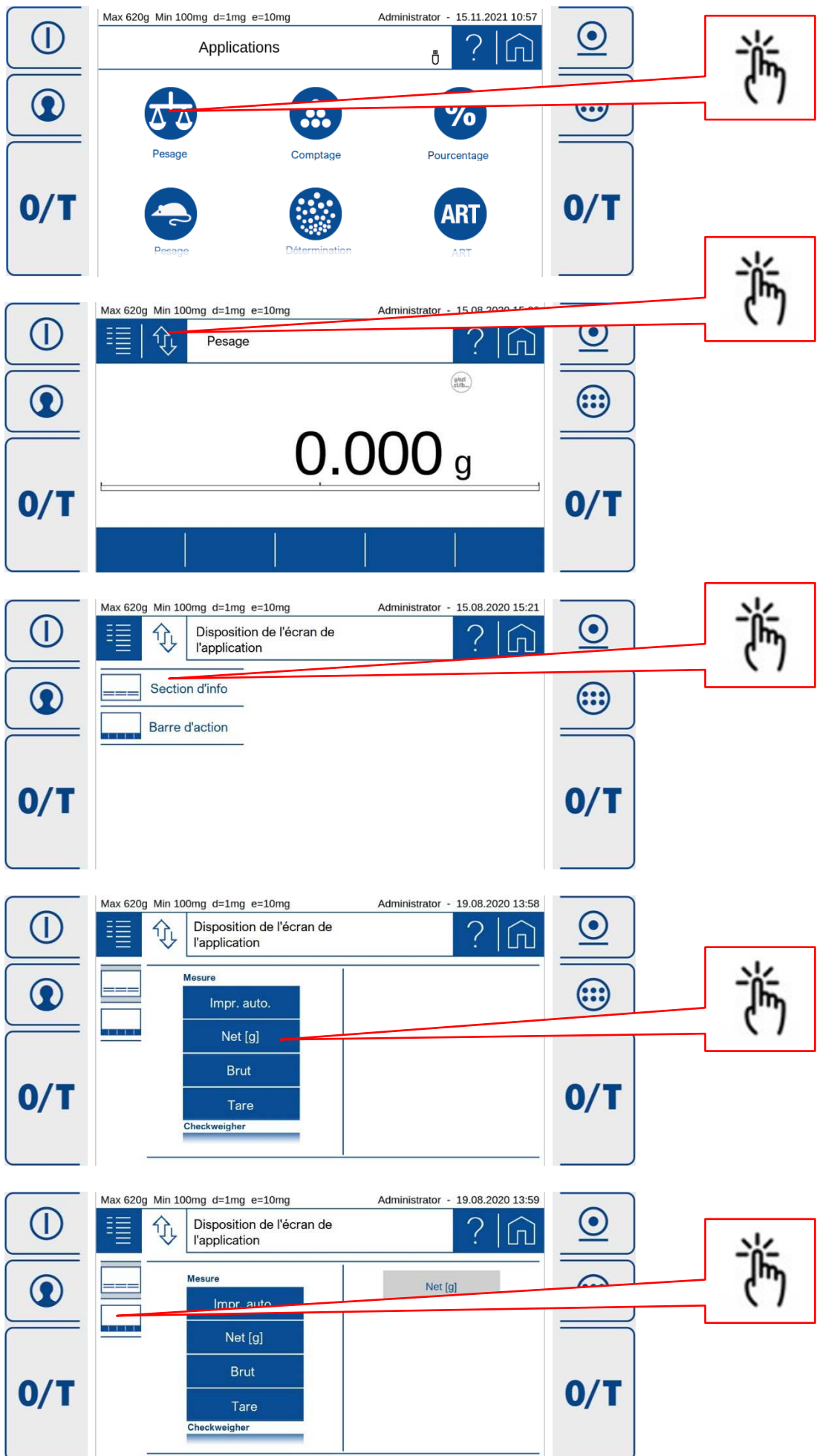
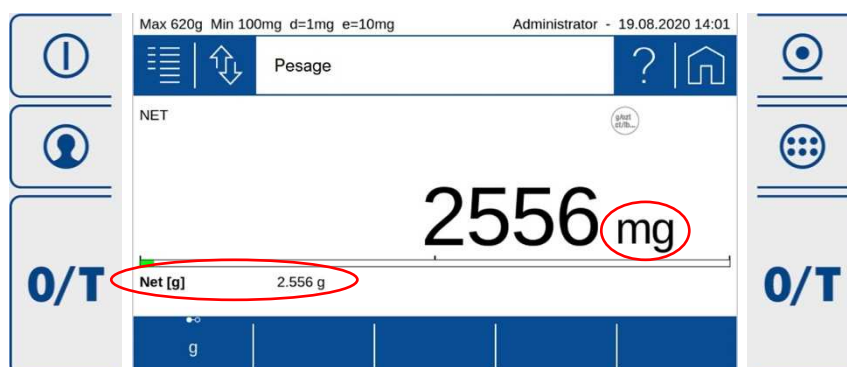
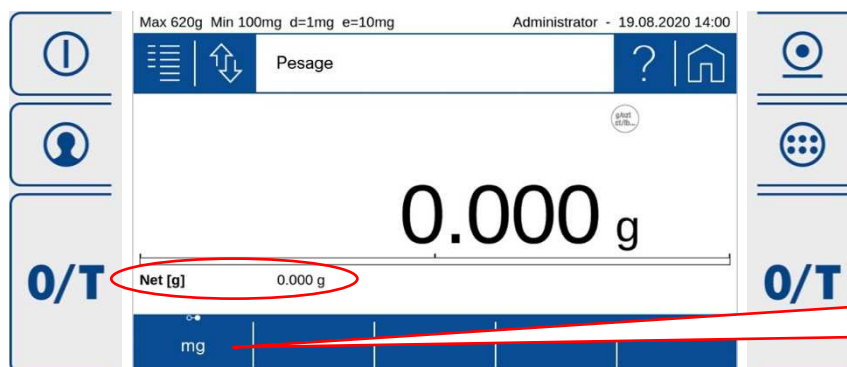
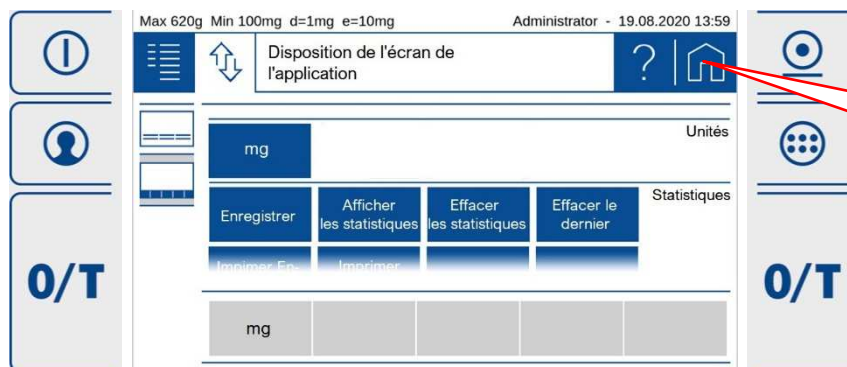
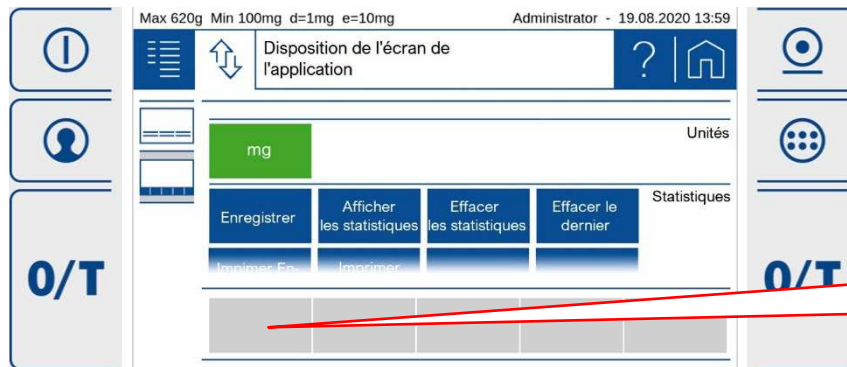
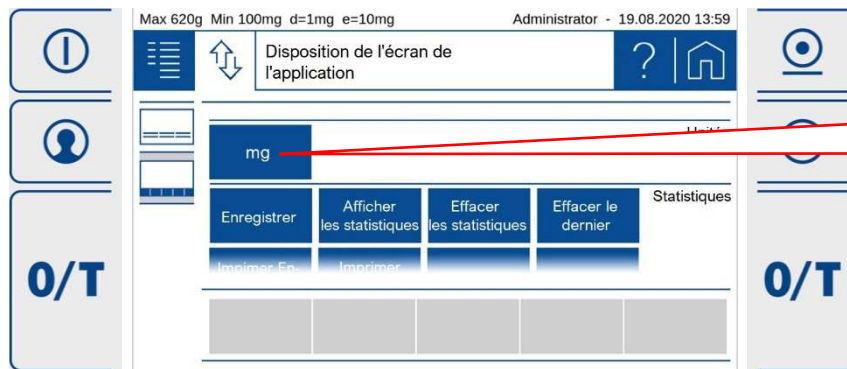
18 VA, 12 V DC, 1.5 A

Note – Hinweis – Remarque – Nota

Retain the original packaging in order to avoid damaging the balance during shipping or transport and to enable the balance to be stored in ideal conditions (5 .. 40°C / 41 .. 104°F) when out of use for long periods.

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No reproduction in any form of this document, in whole or in part, may be made without written authorization from Precisa Gravimetrics AG.

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© Precisa Gravimetrics AG, 8953 Dietikon, Switzerland