

Moisture determination of Rice

This information is intended as a guideline for development of a moisture analysis method and parameter setting.

Materials vary throughout the world and we will be more than happy to test customers' individual samples.



Precisa moisture analysis XM60/120 halogen

Description:	Rice, white powder
Utensils:	Grinder
Preparation:	Ground sample
Sample weight (+/-10%):	2g
Temperature program:	Standard
Temperature:	105°C
Stop-Criteria:	Autostop 2 Digits / 30 s
Moisture content: (average of 6 measurements)	11.79%
Standard Deviation (+/-):	0.03%
Average analysis time:	6.5 min
Reference:	air oven, 105°C for 2 hours 11.88% ± 0.06% (3 measurements)