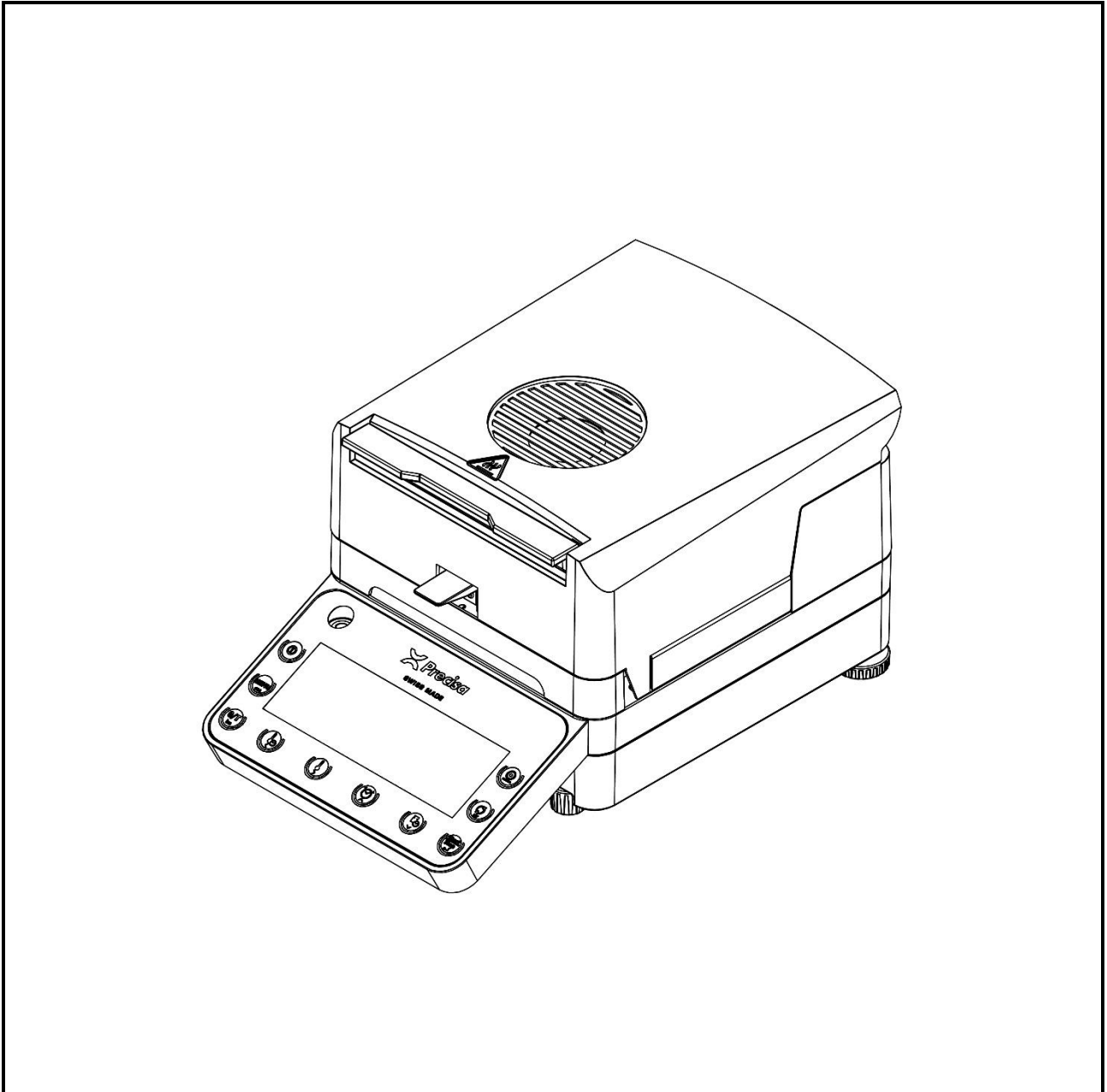
SWISS MADE

# 530 PBM

Quick Start  
Schnelleinstieg  
Départ rapide  
Inicio rápido

M350-8197-900-

 **Precisa**  
THE BALANCE OF QUALITY

## Safety precautions

When the device is used in environments with increased safety requirements, the relevant precautions must be taken.

The device must only be operated with a fault-free power cable.

Before connecting the power cable, ensure that the operating voltage specified on the device matches the mains supply.

The device must only be operated with the power cable supplied.

If safe operation of the device is not possible, it must be disconnected from the power supply and secured against unintentional use.

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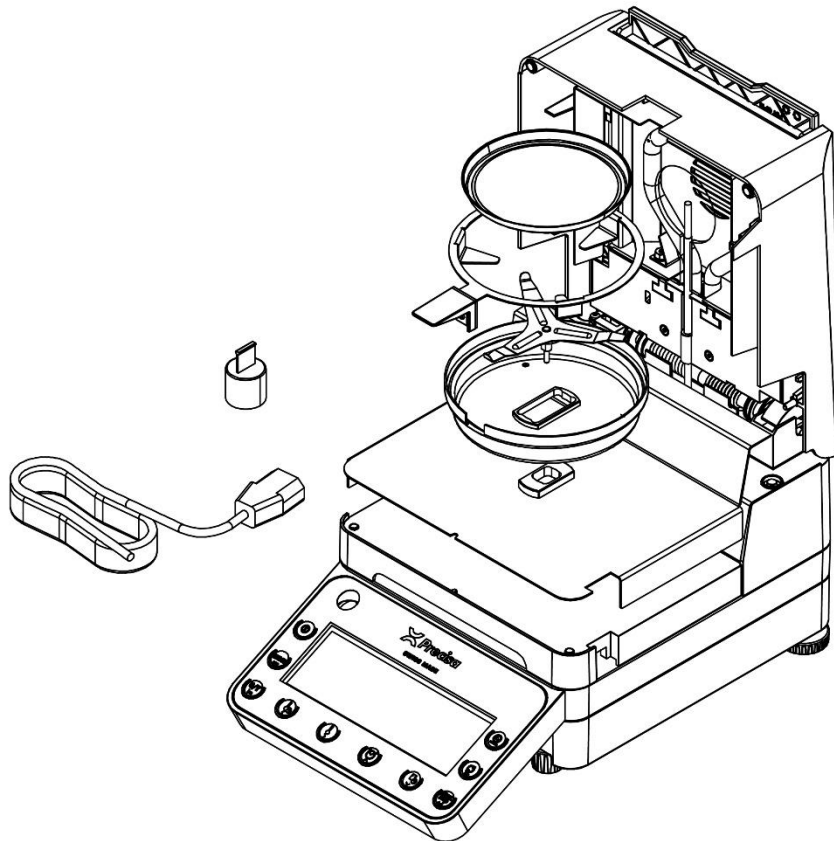
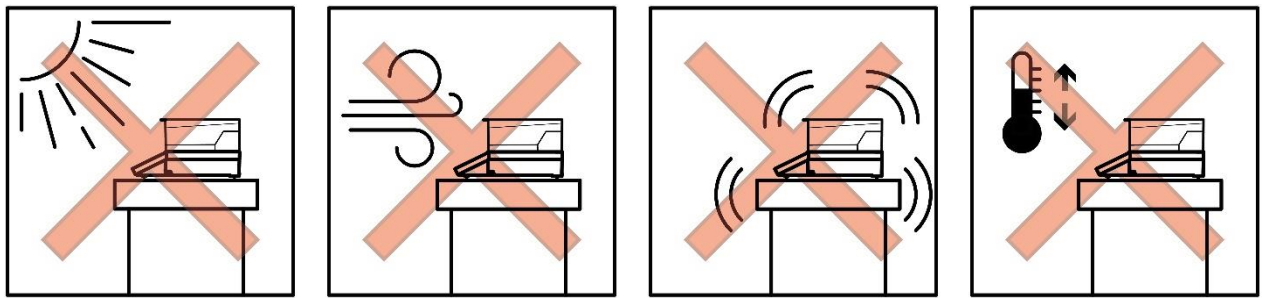
## Starting up – Inbetriebnahme – Mise en service – Puesta en servicio


Only work with a device at operating temperature (warm-up time at least 2 hours). If you have not used the device for a long time, briefly load it with your hand and tare it.

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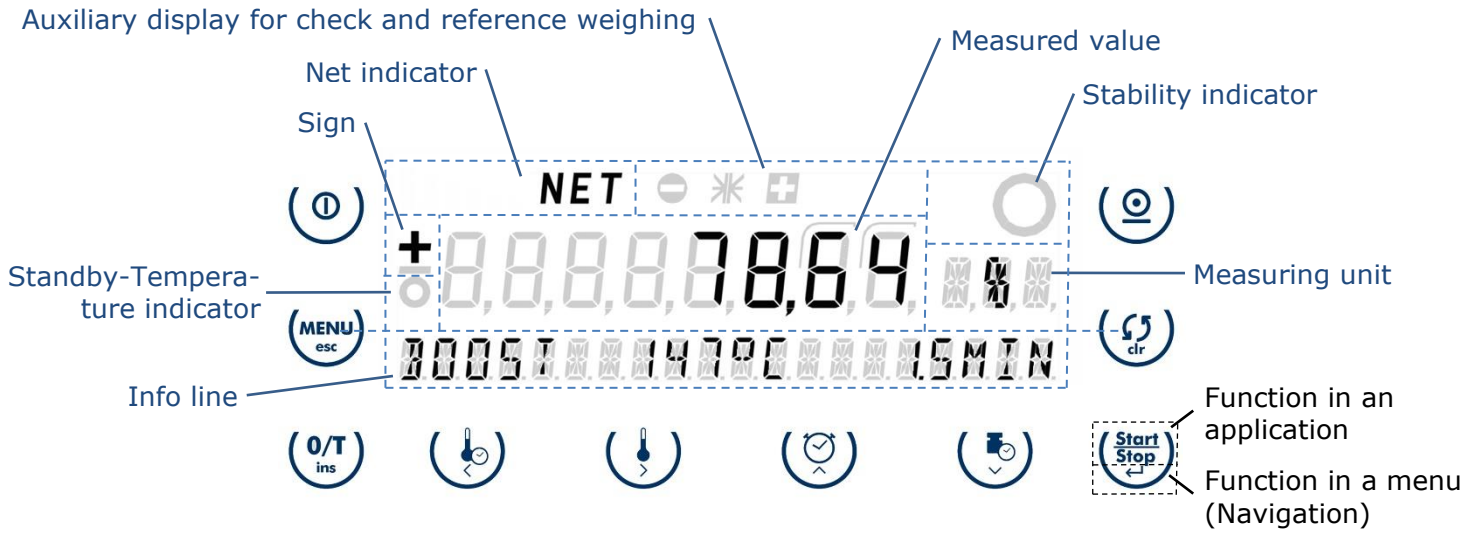[Stop criterium]  
[Down]

[Start/Stop]  
[Enter]

The measured value is stable as soon as all segments of the stability indicator go out. The standby temperature indicator lights up if the standby temperature has not yet been reached.

## The first moisture analysis

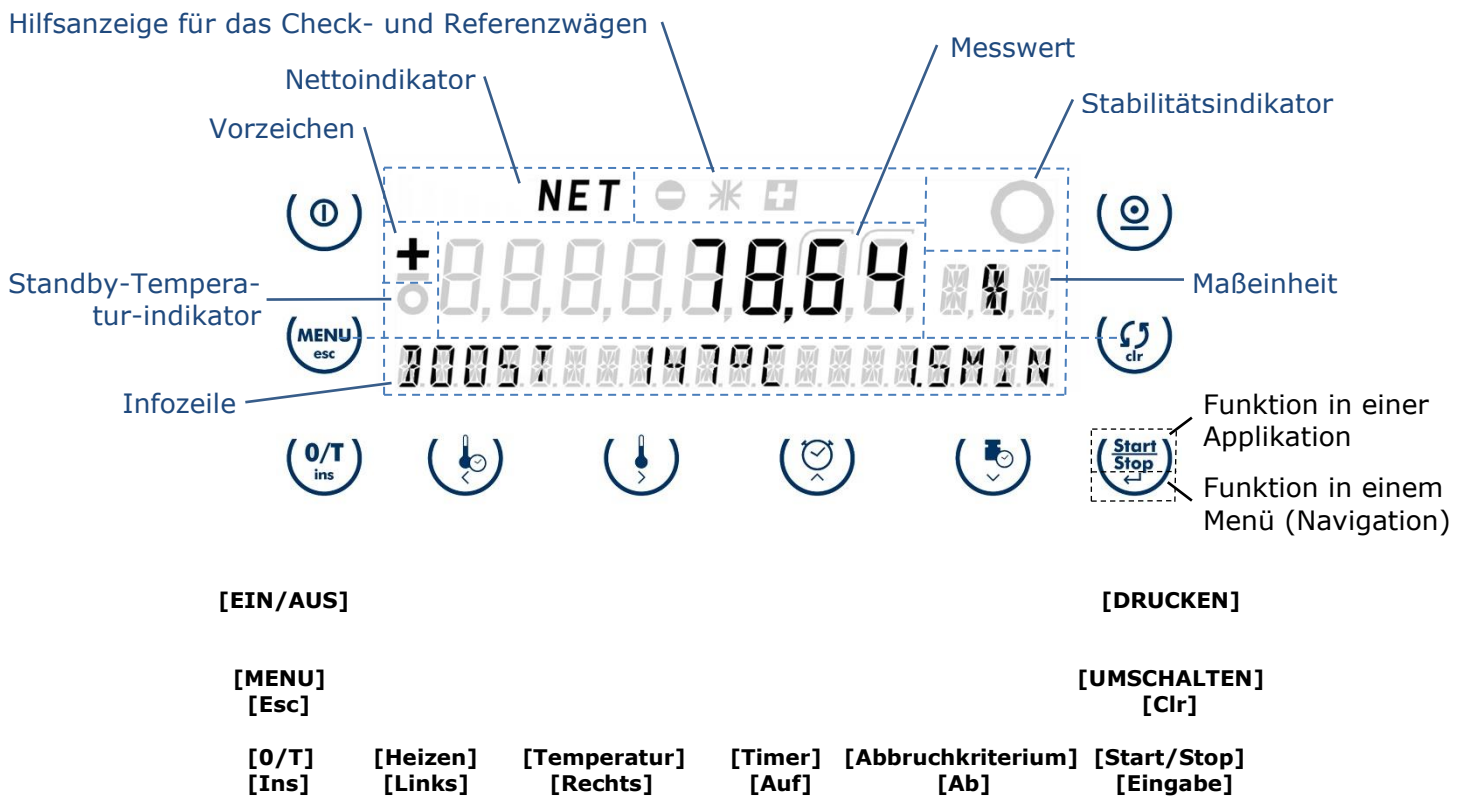
If the device is not set to English: Enter the configuration menu, switch the device off and on with **[ON/OFF]** and, as soon as all segments of the display light up, press and hold **[MODE]** until "SET DATA PRINT" appears. Press **[Down]** repeatedly until "SPRACHE/LANGUE/DIL/IDIOMA" appears. Press **[Enter]**; "SPRACHE/LANGUE/DIL/IDIOMA" starts flashing. Press **[Down]** repeatedly until "LANGUAGE ENGLISH" appears and confirm with **[Enter]**; "LANGUAGE ENGLISH" stops flashing. Press **[Esc]**; the balance finishes starting up.

Press **[Heat]**; "HEAT MODE STANDARD" appears. Press **[Enter]**; "STANDARD" starts flashing. Press **[Down]** repeatedly until "BOOST" appears and confirm with **[Enter]**; "BOOST" stops flashing. Press **[Esc]**. Press **[Timer]**; "STOP TIME 10.0 MIN" appears. Press **[Down]**; "TIMER STOP ON" appears. Press **[Enter]**; "ON" starts flashing. Press **[Down]** repeatedly until "OFF" appears and confirm with **[Enter]**. Press **[Esc]**.

Press **[MENU]** and **[Down]**; "STORE METHOD" appears. Press **[Enter]**; "METHOD \_" appears. Use the cursor keys to enter "DEFAULT\_BOOST" and confirm with **[Enter]**.

Place an empty sample pan and a glass fiber filter on the sample cup support and tare the device with **[O/T]**. Place a paper clip and approx. 1g of water in the sample cup, close the hood and press **[Start/Stop]**; the moisture analysis begins, the display is as in the example above. After a few minutes, the measurement is complete, the hood opens and the result is displayed. Press **[Start/Stop]**; the device is ready for a new measurement.

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© Precisa Gravimetrics AG, 8953 Dietikon, Switzerland.