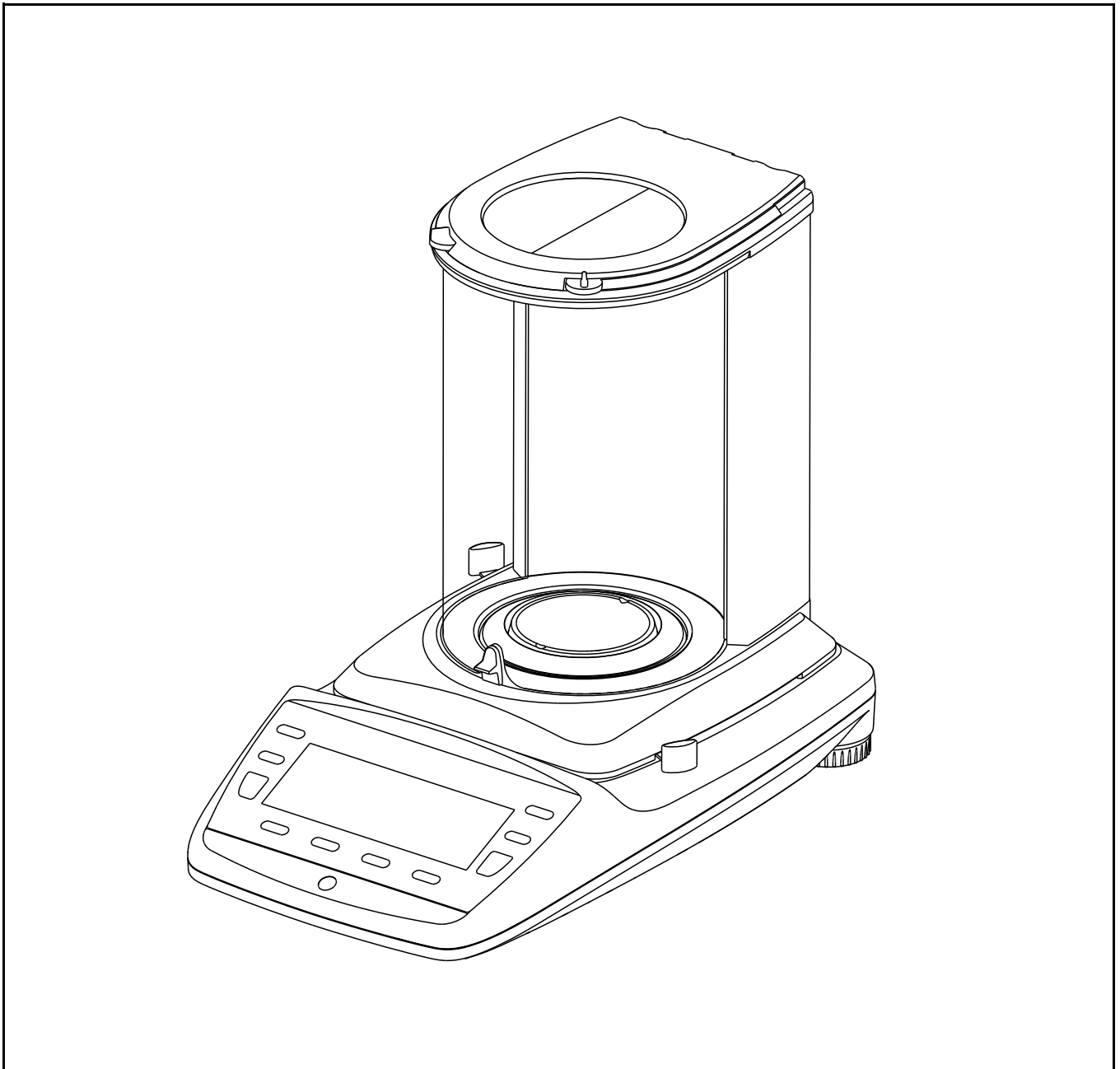
Precisa 360 ES

Brief Operating Instructions
Kurzbedienungsanleitung
Mode d'emploi court

Declaration of conformity

Declaration of conformity for apparatus with CE mark
Konformitätserklärung für Geräte mit CE-Zeichen
Déclaration de conformité pour appareils portant la marque CE
Declaración de conformidad para aparatos con disitintivo CE
Dichiarazione di conformità per apparecchi contrassegnati con la marcatura CE

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Electronic Balance: Precisa Series 360 EP
Precisa Series 360 ES
Precisa Series 321 LS
Precisa Series 321 LX
Precisa Series 321 LT

EU Directive	Standards	EC-type-approval	Issued by
2014/31/EU	EN 45501: 2015	D11-09-010 T7911	PTB NMI

Date: 20.04.2016

Signature:

R. Grolimund R & D Manager

Precisa Gravimetrics AG , Moosmattstrasse 32 , Postfach 352 , CH-8953 Dietikon

1 Safety instructions and Conformity

1.1 Conformity

The device has been manufactured and tested in accordance with the standards and directives set out in the enclosed certificate of conformity.

1.2 Safety instructions

- When using the device in surroundings with increased safety requirements, the corresponding regulations must be observed.
- The device may only be operated with the included power adaptor or power cable.
- Be sure to hold the plug, not the cable, when disconnecting the monitor from an electrical outlet.
- If the power cable, the power adaptor or its cable is damaged, unplug the device immediately from the electrical outlet and keep it from inadvertent operation. The device may only be operated with a power adaptor or a power cable in perfect conditions.
- If there is any reason to believe that it is no longer possible to operate the device without danger, it must be immediately unplugged from the electrical outlet and kept from inadvertent operation.
- In carrying out maintenance work, it is essential to heed the instructions in the chapter "Maintenance and servicing" of the operating instructions.
- Do not operate the device in an area subject to explosion risks or caustic atmospheres.
- Take care when handling with liquids, that no liquid is split into the inside of the balance, on connections of the device, on the power cable or the power adaptor. Should this although happen, unplug the device immediately from the electrical outlet and keep it from inadvertent operation. The device may only be operated again after a re-check made by a service technician.
- The device may only be used for weighing of solid materials, of liquids filled into secure containers, and for running applications designed especially for the device (e.g. animal weighing or density determinations).
- Do not exceed the maximum permissible load of the balance.
- When using the device in combination with other appliances made by the balance manufacturer as well as with appliances produced by other manufacturers, the current regulations for the safe use of the relevant attachments and their application in accordance with instructions must always be observed.

DANGER

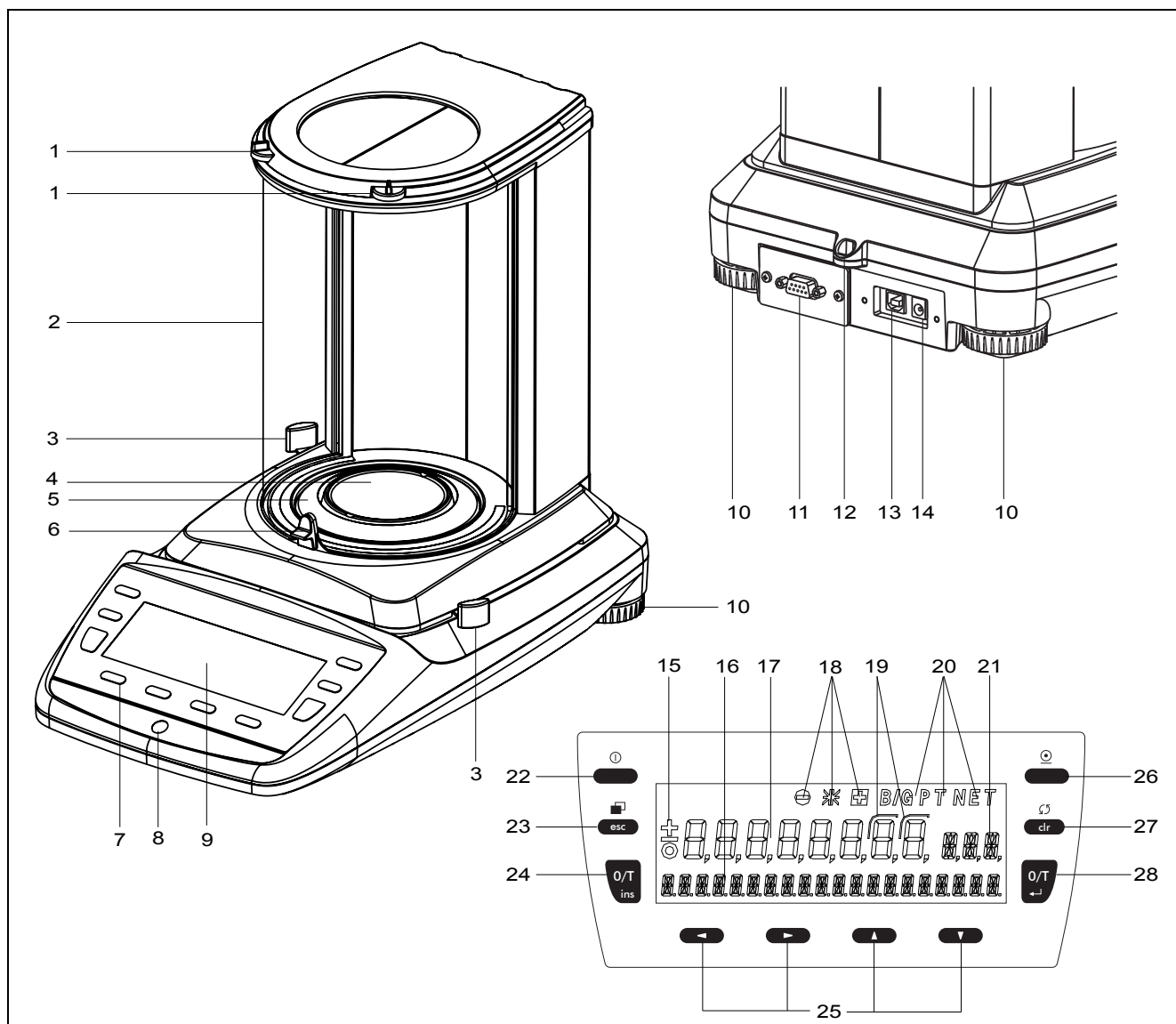
Before connecting the device to the mains, be sure that the operating voltage stated on the device and on the power adaptor or power cable agrees with the local mains voltage. If it doesn't, the equipment must not be connected to the mains at all! In this case contact the Customer Service.

NOTE

If any damage or injury occurs, liability and responsibility rest with the user.

2 Overview

No.	Description	No.	Description
1	Top draft shield opening	15	Prefix display
2	Draft shield	16	Info display
3	Side draft shield opening	17	Measurement display
4	Weighing pan	18	Weight check
5	Protective ring	19	Brackets to mark not approved digits
6	Switch for draft shield opening mode	20	Symbol display
7	10-key control panel	21	Unit display
8	Levelling bubble	22	ON / OFF key
9	Display	23	Menu / ESC key
10	Adjustable feet	24	Tare / Insert key
11	Serial interface, DB9 Female	25	Cursor keys
12	Mechanical anti-theft protection	26	Print key
13	USB device connector for connection to PC	27	Change / Clear key
14	Connecting socket for power adapter	28	Tare / Enter key

3 Operation

3.1 The Keys

3.1.1 Functions of the keys if operated in an application

Key	Designation	Function in Weighing mode
	«ON/OFF»	<ul style="list-style-type: none"> Switching the balance on and off
	«MENU»	<ul style="list-style-type: none"> Calling up the Configuration Menu and the Application Menu
	«T»	<ul style="list-style-type: none"> Initiate Tare Function and/or Calibration Function
	«↻»	<ul style="list-style-type: none"> Switches between the Basic program and the chosen application
	«PRINT»	<ul style="list-style-type: none"> Start print function
	«↑» «←» «→» «↓»	<ul style="list-style-type: none"> Function keys. Start the functions in the info-line.

3.1.2 Functions of the keys if operated in a menu

Keys	Designation	Function in programming mode
	«←», «→»	<ul style="list-style-type: none"> Changes within the menu level
	«↑» «↓»	<ul style="list-style-type: none"> Up/Down movements within the menu Changes selected parameters
	«↵»	<ul style="list-style-type: none"> Selects parameters Stores the changed parameters
	«esc»	<ul style="list-style-type: none"> Interrupts an input Leaves the menu
	«ins»	<ul style="list-style-type: none"> Places insert marker (in text entry)
	«clr»	<ul style="list-style-type: none"> Deletes input (in text entry)
	«PRINT»	<ul style="list-style-type: none"> Inputs a point (in text entry)

3.2 Activating the configuration menu

The basic adjustment of the balance is defined in the configuration.

- Press «**ON/OFF**» to switch the balance on.
- Hold down the «**MENU**» key continuously during the start-up process (which takes about ten seconds) until „UNIT-1“ appears in the display.
- You can now change the configuration menu.

Procedure for changing the language:

Display	Key	Step
	«↓»	Press repeatedly, until the language currently activated is displayed.
	«←»	The language now flashes.
	«↑»	Press repeatedly until the language you require appears.
	«←»	To confirm the choice of language.

This table shows the content of the configuration menu.

Menu	Definable functions
UNIT-1	Unit in which the weighing results are displayed
SET DATA PRINT	Print formats; Type of values to be printed (individual values, continuous printout, time or load change dependent values, date, time, user, etc.)
SET CALIBRATION	Calibration method
SET WEIGHING MODE	Stability mode (Quality of the balance location), Auto-Standby mode, zero correction, tare method (rapid or standard tare)
SET INTERFACE	Baud rate, parity, handshake functions of the peripheral interface
SET DATE AND TIME	Date and time (standard format or American format p.m. and a.m.)
THEFTCODE	Activating/deactivating and changing the anti-theft code.
KEY TONE	Activation of the keypad sound
LANGUAGE	Display language (E, G, F)
CONTRAST	Screen contrast

3.3 Activating the application menu

The balance application programs are called up using the application menu and adapted to the user's needs.

- Press «**MENU**» after the start-up procedure has ended in order to access the application menu.

Menu	Definable functions
SET APP.	Select application program
SETUP APPLICATION	Specify parameters for the application program selected
SET STATISTIC	Statistics and storage functions
SET CHECK +/-	Define nominal weight and limits for comparison weighing
etc.	Other functions available. Refer to the Application Operating Instructions for the description.
AUTO-START ON/OFF	The selected application program can, if required, be loaded automatically every time the balance is switched on
SET USER	Define user profile

■ 3 Operation

3.4 Special operating keys

3.4.1 The tare key

- **Activating taring**
 - Briefly press «T»
- **Range selection**
 - Hold down «T» until „RANGE FINE ON” or „RANGE FINE OFF” is displayed, then release
 - If „RANGE FINE OFF” is selected the balance is working only in the coarse range and therefore the readability is 10 times lower than in the fine range.
- **Activating calibration**
 - Hold down «T» until „CALIBRATION” is displayed, then release
- **Activating an automatic repeatability test (ART)**
 - Hold down «T» until „REPEATABILITY TEST” is displayed, then release

3.4.2 The print key

- **Print out an individual value or a report**
 - Briefly press «PRINT»
- **Reset product counter to 1**
 - Hold down «PRINT» until „RESET PROD.-COUNTER” is displayed, then release
- **Print out a balance status**
 - Hold down «PRINT» until „PRINT STATUS” is displayed, then release
- **Print out the application-setup**
 - Hold down «PRINT» until „PRINT APPLICATIONS” is displayed, then release
- **Print out the calibration information**
 - Hold down «PRINT» until „PRINT CALIBRATIONS” is displayed, then release «PRINT»

3.4.3 The change key

- **Switch to other applications**
 - As long as you hold «», all active applications are shown one after the other: if, for example, the statistics program, the check program and the count application are activated, „WEIGHING”, „STATISTIC”, „CHECK” and „COUNT” appear in the info-line one after the other:
 - Release «» when the application to which the balance should be switched, appears in the display.

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