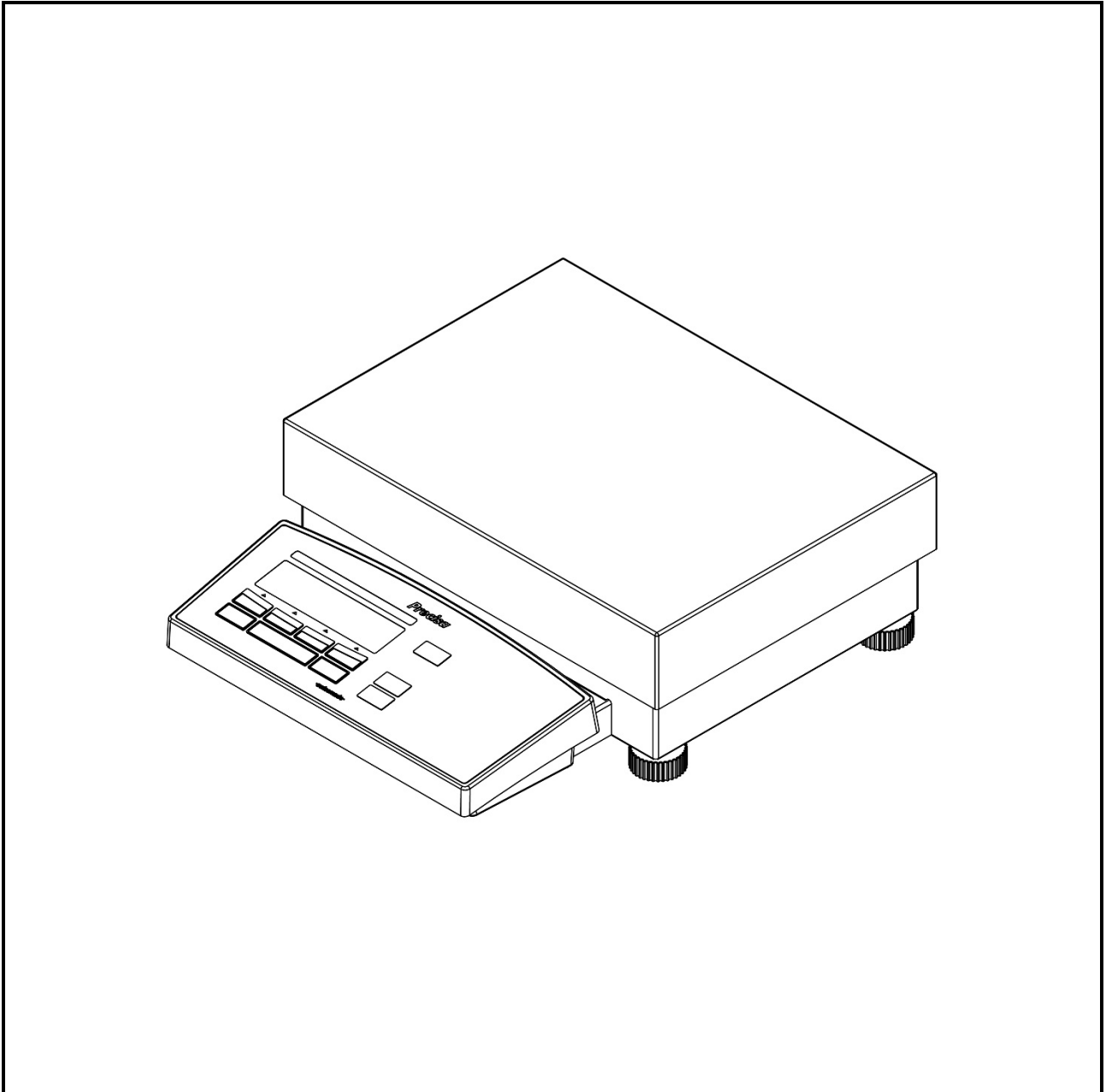
SWISS MADE

# 490 IBK/ITK/ISK

Quick Start  
Schnelleinstieg  
Départ rapide  
Inicio rápido

M350-8191-900-

 **Precisa**  
THE BALANCE OF QUALITY

## Safety precautions

When the balance is used in environments with increased safety requirements, the relevant precautions must be taken.

The balance must only be operated with a fault-free power adapter.

Before connecting the power adapter, ensure that the operating voltage specified on the power adapter matches the mains supply.

The balance must only be operated with the power adapter supplied.

If safe operation of the balance is not possible, it must be disconnected from the power supply and secured against unintentional use.

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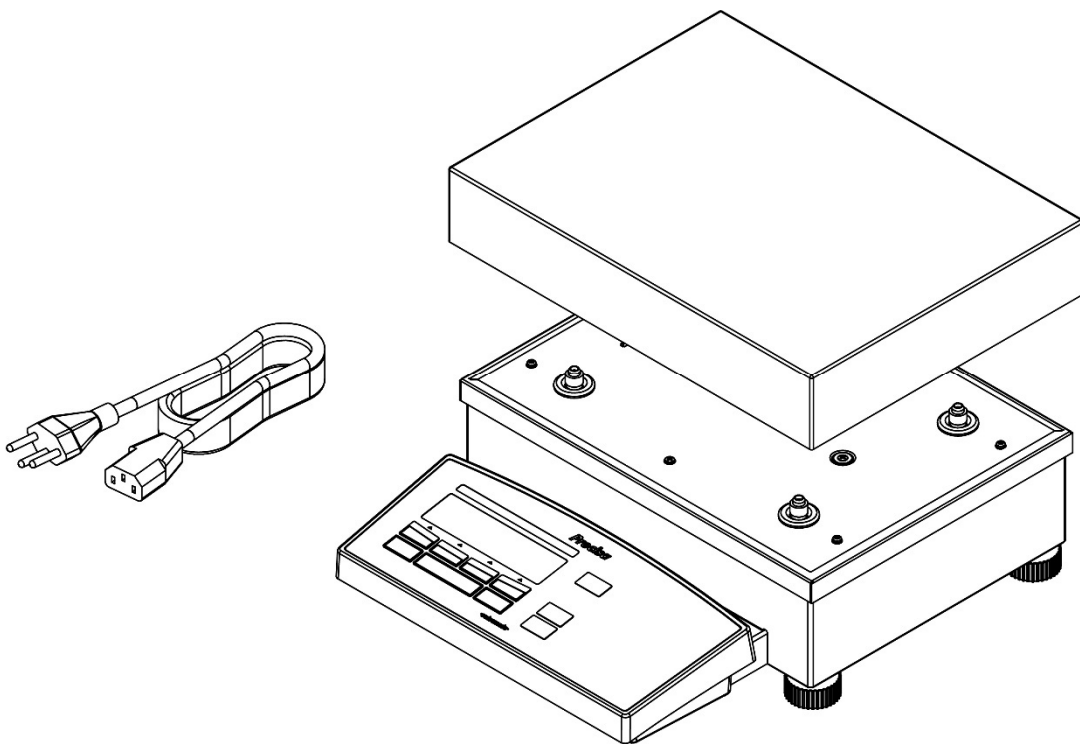
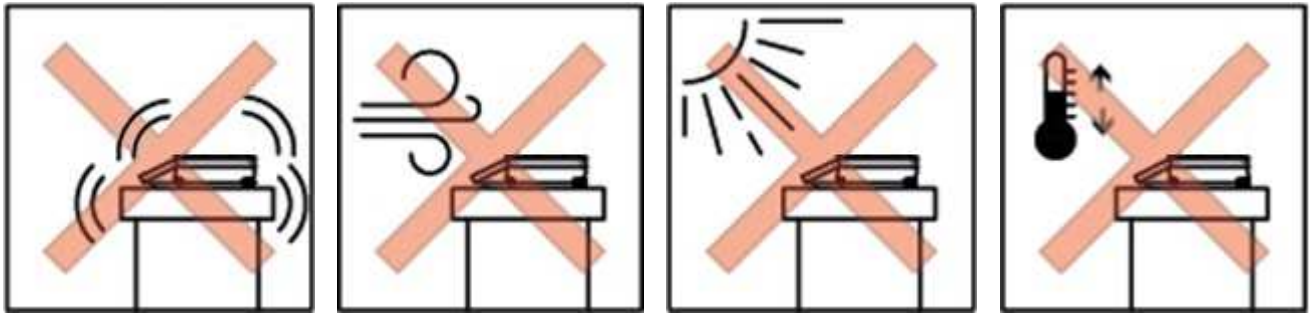
## Starting up – Inbetriebnahme – Mise en service – Puesta en servicio


Only work with a balance at operating temperature (warm-up time at least 2 hours). If you have not used the balance for a long time, briefly load it with your hand and tare it.

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## Operation

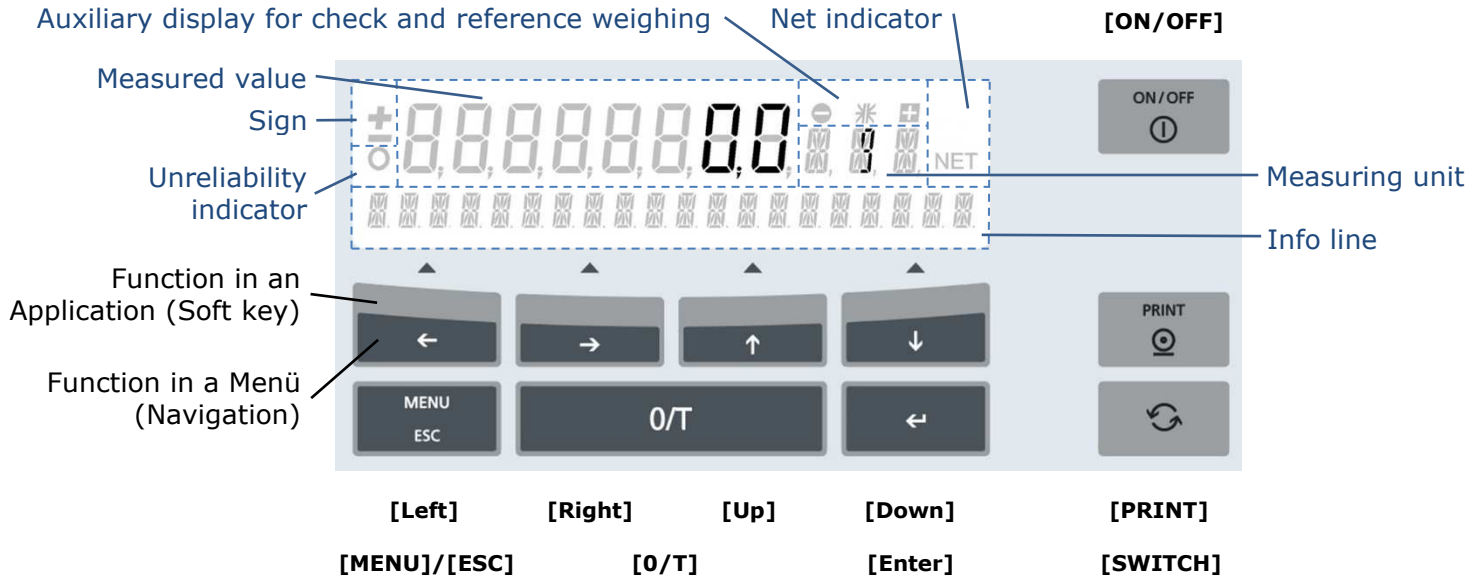The measured value is stable as soon as the unit, e.g. "g", appears. The unreliability indicator lights up while the scale is recalculating the reference weight, for example, during counting, or when it is verified and in the warm-up phase.

## Application example: Adding up components (application net total)

If the scale is not set to English, enter the configuration menu: Switch the scale off and on with **[ON/OFF]** and, as soon as all segments of the display light up, press and hold **[MENU]** until "UNIT-1 g" appears. Press **[Down]** repeatedly until "SPRACHE DEUTSCH/LANGUE FRANÇAISE" appears. Press **[Enter]**; "SPRACHE DEUTSCH/LANGUE FRANÇAISE" starts flashing. Press **[Down]** repeatedly until "LANGUAGE ENGLISH" appears and confirm with **[Enter]**; "LANGUAGE ENGLISH" stops flashing. Press **[ESC]**; the scale finishes starting up.

To enter the application menu, press **[MENU]**; "SET APP. OFF" appears. Press **[Enter]**; "OFF" starts flashing. Press **[Down]** repeatedly until "NET-TOTAL" appears and confirm with **[Enter]**; "NET-TOTAL" stops flashing. Press **[ESC]**; the scale exits the application menu, starts the application Net total and displays 0g as the measured value and "STO 0 RES INF" in the info line.

Place the first component, e.g. 1000g, and press **{STO}**; the scale records the value, tares itself and displays "STO 1 RES INF" in the info line. Place the second component, e.g. 2000g, and press **{STO}**, and finally place the third component, e.g. 200g, and press **{STO}**.

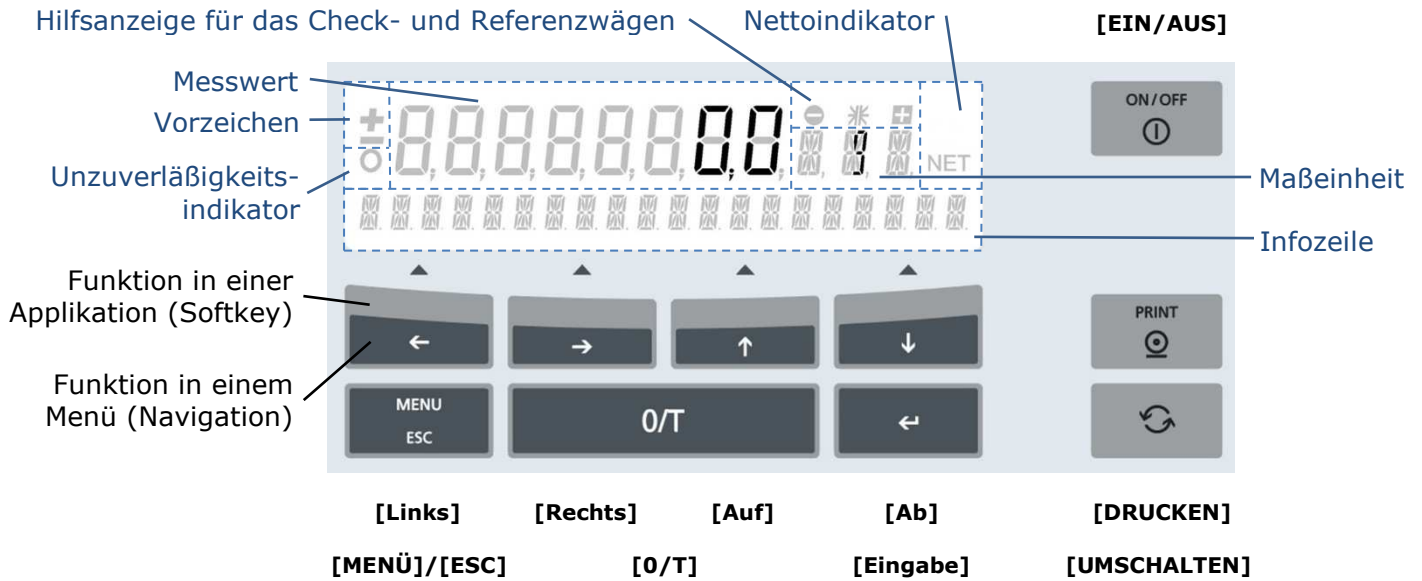
While adding up, the current total weight can be viewed at any time by pressing **[SWITCH]**. The scale then switches to weighing mode. To return to the application, press **[SWITCH]** again. To switch from the analogue display to the application commands in the info line, press **[SWITCH]** twice; the scale switches to the weighing mode and back to the application. It then displays e.g. "STO 3 RES INF".

Press **{INF}**; the scale displays in the info line all information on the weights just added up. Use **[Down]** and **[Up]** to scroll through the entire stack of info lines: Total, remaining capacity and all individual weights, e.g. 1000g, 2000g and 200g. Press **[ESC]** to exit the information query; the scale switches back to the running application and is ready for another weighing-in.

To restart, press **{RES}**; the scale resets the counter to 0.

To quit the application, press **[SWITCH]**; the scale switches back to weighing mode.

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© Precisa Gravimetrics AG, 8953 Dietikon, Switzerland.