

Release Notes 1.5.0

New features

- Alibi storage available for application Weighing

Miscellaneous

- EDP mode, revised

Bug fixes

- Audit trail, date search
- Identification, show/hide soft button
- Adjustment history, deviation calculation

Release Notes 1.4.0

New features

- First **EU/OIML/UKCA** approved version. (T12340)

Miscellaneous

- Manual adjustment, permitted for Standard and Restricted users.

Bug fixes

- Remote network communication corrected.

Release Notes 1.3.0

New features

- Add-on Peak-Hold available for application Weighing
- Counting, reference optimization added

Miscellaneous

- Protocol, print item "Form feed" added
- Protocol, HEX character entry for "Text paragraph" added
- Counting, soft button for piece weight entry introduced
- Counting, soft button for individual piece entry introduced
- Language Polish added

Bug fixes

- HID, wrong date string corrected
- Frozen screen when displaying message boxes

Release Notes 1.2.1

Bug fixes

- Minimum weight indicator (C, D and G balances)
- Zero setting range (approved balances)

Release Notes 1.2.0

New features

- Density application
- Add-on Checkweigher available for application Weighing, Counting, Percent and Dynamic weighing
- Wifi connection by USB Wifi adapter (350-8941)

Miscellaneous

- Soft button for levelling added
- Soft button for preset tare added
- Weight entries for the current application in the selected application unit
- Reference unit added to the System Settings menu for global weight entries
- Unit kilogram and pound added
- Info item Gross and Tare added
- Item for protocol format length 120 added
- Returning from Adjustment or Levelling to the main application by [Home] button
- Linearization with displayed default values

Bug fixes

- Unit Bath conversion factor corrected
- Unit Tola conversion factor corrected
- Unit Teal Singapore corrected
- Dynamic weighing tare/zero function corrected