

Release Notes 1.3.5

New features

- Context help in French language included

Miscellaneous

Bug fixes

- Possible blockage when turning on fixed (issue in 1.3.2, 1.3.3)
- Minor bug fixes

Release Notes 1.3.4

New features

- Protocol type "SICS V1.0" included.
Remote command list included in context help of setting "Protocol" and "Data exchange - ..."
- Checkweigher: Signal lamp control on RS232 serial port included.

Miscellaneous

- Max. number of users increased to 25

Bug fixes

- Possible blockage when turning on fixed (issue in 1.3.2, 1.3.3)
- Translated texts corrected or optimized
- Minor bug fixes

Release Notes 1.3.3

New features

Miscellaneous

- Remote command list included in context help of setting "Data exchange - ..."
- Soft button "1/...d" renamed to "x Decimals"

Bug fixes

- Info item "ID": Text-initialisation when changing application

Release Notes 1.3.2

New features

- Application "Dynamic" add-on "Checkweigher" included
- Soft button "1/...d" for resolution-reduction included

Miscellaneous

- Add-on "Peak Hold" protocol item "Peak value" included
- Add-on "Statistics" action button "Max. n" included
- MSW setting "Capture below minimum" included
- For Statistics- and Protocol-Load change adjustable "Min. Load" included
- Data exchange mode "Measurement transfer" transfer item "ID" included
- Remote control commands "D<text>", "DN" and "@<text>", "@N" for writing text into the display included
(Remote command list is now included in context help "Data exchange – Serial Port")

Bug fixes

- French text corrected

Release Notes 1.3.1

New features

- Printer LP4024, printout in Chinese language available (no settings on printer required).
Note: Printer Firmware LP 4024 Vers:3.0 required

Max 225g Min 1mg d=0.01/0.1mg e=1mg Administrator - 22.05.2024 16:02

系统设置		?	🏠
称重参数	模式	协议	
数据交换 ▲	波特率	9600	
串口	奇偶校验	8-无-1	
协议 ▼	握手协议	无	
重量转换	格式化	字符串40	
外设设置	打印机	LP4024	

通用设置

Miscellaneous

- Translations revised
- Context help for addon Identification added

Bug fixes

- Add-On Identification: Customizing of soft buttons [ID1..6]
- Add-On Checkweigher: Function of soft button [Capture nominal]

Release Notes 1.3.0

New features

- Time synchronisation via the network added, works from terminal version BSP V2.2.4 upwards (see device information).
- Time zone setting added. Automatic daylight-saving time regulation included.

Miscellaneous

- Data output (serial port, mass storage and network) in Chinese language encoded as character set GB18030 for traditional Chinese and GB2312 (subset of GB18030) for simplified Chinese.

Bug fixes

Release Notes 1.2.4

New features

Miscellaneous

Bug fixes

- RTC deviation correction

Release Notes 1.2.3

New features

Miscellaneous

- Service remote command CLRHISTORYcccc added.
- Public remote command FWHISPRINT added.

Bug fixes

Release Notes 1.2.1

New features

Miscellaneous

- Writing to a CSV file revised. CSV separator selectable, semicolon or comma

Bug fixes